INSTRUCTION MANUAL

Analytical Balance

GH-120

GH-200

GH-300

GH-202

GH-252



© 2008 A&D Company Ltd. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company Ltd.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.

Windows, Word and Excel are registered trademarks of the Microsoft Corporation.

Contents

Basic Op	peration	
1. 1.1. 1.2. 1.3.	Introduction About This Manual Features Compliance	4
2. 2.1.	Unpaking the BalanceInstalling the Balance	
3.1. 3.2. 3.3. 3.4.	Precautions Before Use During Use After Use Power Supply	8
4.	Display Symbols and Key Operation	10
5. 5.1. 5.2.	Weighing UnitsUnitsChanging the Units	11
6. 6.1. 6.2. 6.3. 6.4.	Weighing Basic Operation (Gram Mode) Dual Range Counting Mode (PCS) Percent Mode (%)	16 17 18
Adaptin	ng To The Environment	
7. 7.1. 7.2.	Response Adjustment	21
8. 8.1. 8.2. 8.3. 8.4. 8.5. 8.6. 8.7.	Calibration Calibration Group Automatic Self Calibration One-Touch Calibration Calibration Test Using the Internal Mass Calibration Using an External Weight Calibration Test Using an External Weight Correcting the Internal Mass Value	
9. 9.1. 9.2.	Function Switch and Initialization Permit or InhibitInitializing the Balance	29
Selectin	g Functions	
10. 10.1. 10.2. 10.3. 10.4.	Function Table	34 34 37
10.5.	Description of the Item "Data Format"	

10.6.	Data Format Examples	
10.7.	Clock and Calendar Function	44
11.	ID Number and GLP Report	46
11.1.	Setting the ID Number	46
11.2.	GLP Report	47
12.	Data Memory	52
12.1.	Notes on Using Data Memory	52
12.2.	Data Memory for Weighing Data	
12.3.	Data Memory for Calibration and Calibration Test	56
13.	Underhook	57
14.	Density Measurement	58
Interface	e And Communication	
15.	Standard Input and Output Interface	61
15.1.	RS-232C Interface	
15.2.	Connection to Peripheral Equipment	
15.3.	Commands	
Mainten	ance	
16.	Maintenance	70
16.1.	Treatment of The Balance	
16.2.	Error Codes	
16.3.	Other Display	
16.4.	Checking the Balance Performance and Environment	
16.5.	Asking for Repair	
17.	Specifications	
17.1.	External Dimensions	
17.2.	Options and Peripheral Equipment	76
18.	Terms/Index	80
18.1.	Terms	80
18.2.	Index	81

1. Introduction

This manual describes how the balances of GH series work and how to get the most out of them in terms of performance. Read this manual thoroughly before using the balance and keep it at hand for future reference.

1.1. About This Manual

1.2. Features

- Automatic self calibration, using the internal mass, adapting to changes in temperature.
- Response adjustment adapting to drafts and/or vibration automatically.
- Memory function to store weighing data and calibration data.
 When weighing data is stored only, 200 data can be stored in maximum.
 Interval memory mode to store the weighing data periodically.
- Good laboratory practice (GLP) data can output using the RS-232C serial interface.
- A built-in clock and calendar that can add the time and date to the output data.
- Underhook, for measuring density and weighing magnetic materials.
- Multiple weighing units with most of the common units used around the world.
- Density mode for calculating the density of a solid.
- Standard RS-232C serial interface to communicate with a computer.
 Windows communication tools software (WinCT) to allow easy communication with a computer using Windows.
- When multiple balances have the GH-08 Ethernet interface installed and are connected to a LAN, data can be acquired from them using the WinCT-Plus software.

1.3. Compliance

1.3.1. Compliance With FCC Rules

Please note that this equipment generates, uses and can radiate radio frequency energy. This equipment has been tested and has been found to comply with the limits of a Class A computing device pursuant to Subpart J of Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when equipment is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference.

(FCC = Federal Communications Commission in the U.S.A.)

1.3.2. Compliance With EMC Directives

This device features radio interference suppression and safety regulation in compliance with the following Council Directives

Council directive 89/336/EEC EN61326 EMC directive

Council directive 73/23/EEC EN60950 Safety of Information Technology Equipment

The CE mark is an official mandatory European marking.
 Please note that any electronic product must comply with local laws and regulations when sold or used anywhere outside Europe.



CE

A & D Instruments Ltd. hereby declare that the following weighing products conform to the requirements of the council directives on ...

Electromagnetic Compatibility (EMC) 89/336/EEC and Low Voltage Equipment (LVD) 73/23/EEC amended by 93/68/EEC

provided that they bear the CE mark of conformity as shown above.

GH and HR-i Series

Standards applicable:

EN 61326 1997 +A1: 1998, +A2: 2001 Class B Electrical equipment for measurement, control and

laboratory use - EMC requirements for emission and

minimum requirements for immunity.

Specification for safety of household and similar electrical appliances. LVD General requirements EN 60335-1:1991

Safety of Information Technology Equipment. EN 60950

CE Mark first applied October 2004

Warning

These may be class A products. In a domestic environment these products may cause radio interference in which case the user may be required to take adequate measures.

Signed for A&D Instruments in Oxford England October 2005

Takeo Goto Managing Director








2. Unpacking the Balance


Unpack the balance carefully. Keep the packing material to be used for transporting the balance in the future. See the illustrations to confirm that everything is contained.

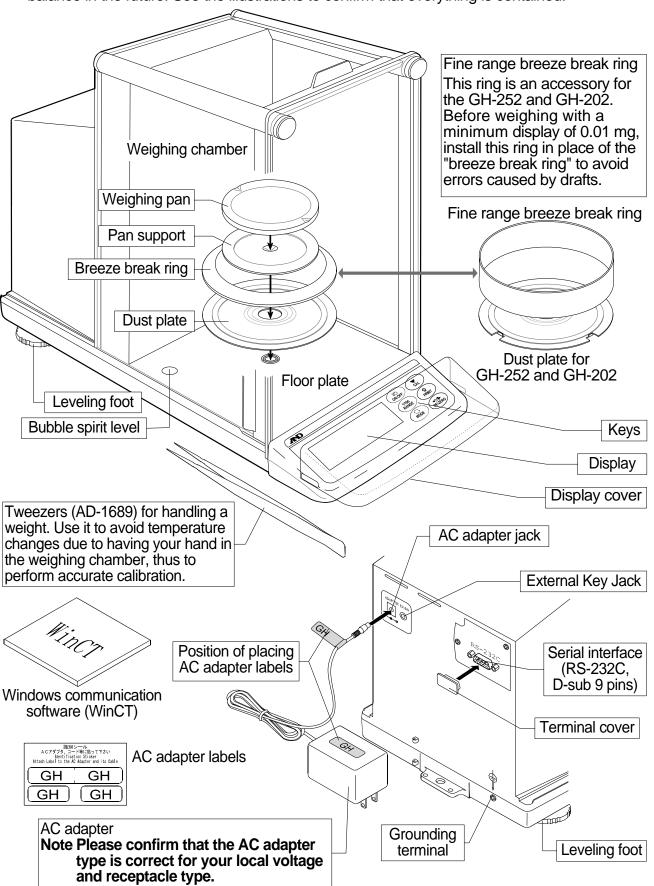


2.1. Installing the Balance

Install the balance as follows:

- 1 Consider the section "3. Precautions " for installing your balance. Place the balance on a firm weighing table.
- 2 Assemble the "Dust Plate", "Breeze Break Ring" (or "Fine range breeze break ring") "Pan support", and "Weighing Pan" on your balance. There is a reference illustration on the previous page.
- 3 Adjust the level of the balance using the leveling feet. Ground the balance chassis for discharging static electricity if you have a static problem.
- 4 Please confirm that the adapter type is correct for your local voltage and power receptacle type.
- 5 Connect the AC adapter to the balance. Warm up the balance for at least one hour with nothing on the weighing pan.
- 6 Calibrate the balance before use. Refer to "8. Calibration".

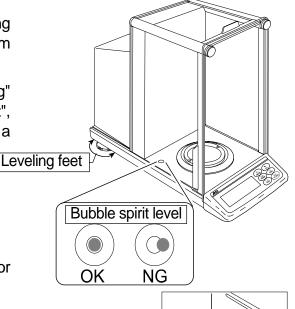

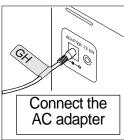

3. Precautions

3.1. Before Use

To ensure that you get the most from your balance, please try to follow these conditions as closely as possible.


- Please confirm that the AC adapter type is correct for your local voltage and receptacle type.
- Ensure a stable power source when using the AC adapter.
- □ The best operating temperature is about 20°C / 68°F at about 50% Relative Humidity.
- The weighing room should be free of dust.
- The weighing table should be solid and free from vibration, drafts (such as frequently opening doors or windows) and as level as possible.
- Corners of rooms are best as they are less prone to vibration.
- Do not install the balance near a heater or air conditioner.
- Do not install the balance in direct sunlight.
- Do not use the balance near other equipment which produces magnetic fields.
- Adjust the level of the balance using the leveling feet.
- Please warm-up the balance for at least one hour. Plug-in the AC adapter as usual.
- Calibrate the balance before using and after moving it to another location.


 \triangle Do not place or use the balance where there is flammable or corrosive gas present.

3.2. During Use

Note the following items to get accurate weighing data.

- Discharge static electricity from the weighing material.
 When weighing sample (plastics, insulator, etc.) could have a static charge, the weight value is influenced.
 Ground the balance, and
 - Eliminate the static electricity using the optional AD-1683.
 - Try to keep the ambient humidity above 45%RH at the room.
 - Use a metal shield case.
 - Wipe a charged material (plastic sample etc.) with a damp cloth.

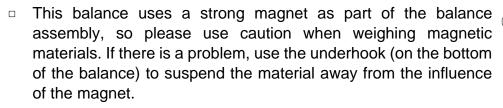

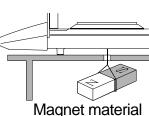


Metal case

Grounding

Charged material

Eliminate temperature differences between the weighed sample and the environment. When a sample is warmer (cooler) than the ambient temperature, the sample will appear lighter (heavier) than the true mass. This error is due to the rising (falling) draft next the sample. If you touch the sample, the same type error will occur. Do not touch the sample directly with your hand. Use tweezers or other tools.

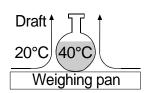

- Make each weighing gently and quickly to avoid errors due to changes in the environmental conditions.
- Before weighing with a minimum display of 0.01 mg for the GH-252 and GH-202, the "fine range breeze break ring" can be installed in place of the "breeze break ring" to avoid errors caused by drafts.



 Do not drop things upon the weighing pan, or place a weight beyond the range of the balance on the weighing pan.



Do not use a sharp instrument (such as a pencil or ball point pen) to press the keys, use your finger only.






- Press the RE-ZERO key before each weighing to prevent possible errors.
- Calibrate the balance periodically so as to eliminate possible errors.
- Take into consideration the affect of air buoyancy on a sample when more accuracy is required.
- Avoid foreign matter (dust, liquid or metal fragments) that could get inside the balance.
- Operate your balance gently. Shorten the operation time as much as possible (Opening and closing door, putting and removing sample). Use a pair of tweezers (pincette) to avoid temperature changes due to heat from inserting your hand into the weighing chamber.



3.3. After Use


- Avoid mechanical shock to your balance.
- Do not disassemble the balance. Contact your local A&D dealer if your balance needs service or repair.
- Do not use solvents to clean the balance. For best cleaning, wipe with a dry lint free cloth or a lint free cloth that is moistened with warm water and a mild detergent.
- Avoid foreign matter (dust, liquid or metal fragments) that could get inside the balance.

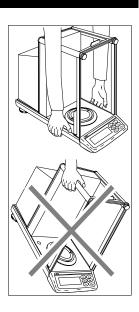

3.4. Power Supply

- Do not remove the AC adapter while the internal mass is in motion, for example, right after the AC adapter is connected, or during calibration using the internal mass. If the AC adapter is removed under the conditions described above, the internal mass will be left unsecured, that may cause mechanical damage when the balance is moved. Before removing the AC adapter, press the ON:OFF key and confirm that zero is displayed in weighing mode.
- When the AC adapter is connected, the balance is in the standby mode if the standby indicator is on. This is a normal state and does not harm the balance. For accurate weighing, we recommend that you always plug in your balance so it can warm up.

4. Display Symbols and Key Operation

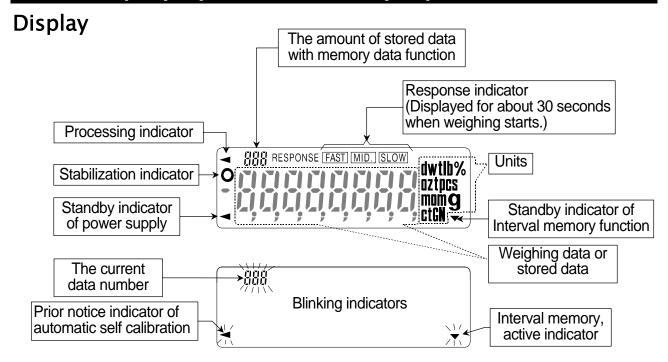

Key operation

Press and release the key immediately" or "Click the key"



□ Press and hold the key ······




Key	When pressed and released	When pressed and held					
I/O ON:OFF	The key to turn the display ON and OFF. The standby indicator is displayed when the display is turned off. The weighing mode is enabled when the display is turned on. This key is available anytime. Pressing the key during operation will interrupt the operation and turn the display OFF.						
1/10d RANGE	 In the weighing mode, the key to turn the minimum weighing value ON and OFF. In the counting or percent mode, the key to enter the sampling mode. 	The key to enter the function table mode. Refer to "10. Function Table".					
MODE	The key to switch the preset weighing units stored in the function table. Refer to "5. Weighing Units".	The key to perform automatic response adjustment.					
CAL	The key to perform calibration using the internal mass.	The key to display other items of the calibration menu.					
Q PRINT	The key to outputs the weighing data to a printer or personal computer (or store it in memory) depending on the function table settings. (Factory setting = output)	No function. (factory setting) By changing the function table: "Title block" and "End block" for GLP report are output. The data memory menu is displayed.					
→0/T+ RE-ZERO	The key to set the display to zero.						

5. Weighing Units

5.1. Units

All weighing units and weighing modes are as follows:

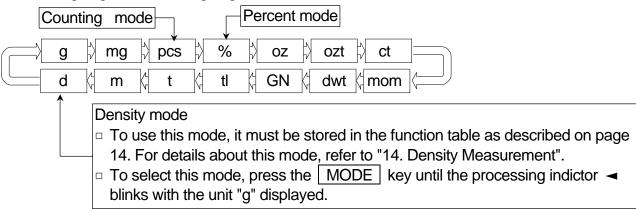

- A unit or mode can be selected and stored in the function table as described in "5.2.
 Changing the Units".
- If the law in your area permits, you may use all of the units. You can disable the units that you don't regularly use. And you are able to turn them back on.
- If a weighing mode (or unit of mass) has been turned off, that mode or unit will be missing in the sequence. Tael has four varieties, one of which can be selected and installed at the factory.
- □ To select a unit or mode for weighing, press the MODE key.
- For details about the units and modes, see the table below:

Name (unit, mode)	Abbreviation	Display unit	Conversion factor	
Gram	g	g	1 g	
Milligram	mg	mg	0.001 g	
Counting mode	pcs	pcs	-	
Percent mode	%	%	-	
Ounce (Avoir)	OZ	OZ	28.349523125 g	
Troy Ounce	ozt	ozt	31.1034768 g	
Metric Carat	ct	ct	0.2 g	
Momme	mom	mom	3.75 g	
Pennyweight	dwt	dwt	1.55517384 g	
Grain (UK)	GN	GN	0.06479891 g	
Tael (HK general, Singapore)			37.7994 g	
Tael (HK jewelry)	tl	tı	37.429 g	
Tael (Taiwan)	u	LI LI	37.5 g	
Tael (China)			31.25 g	
Tola (India)	t	t	11.6638038 g	
Messghal	m	m	4.6875 g	
Density mode	d	d Refer to "14. Density Measurement"		

The tables below indicate the weighing capacity and the minimum display for each unit, depending on the balance model.

	GH-252						
Unit	Standar	d range	Precision range				
Offic	Capacity	Minimum display	Capacity	Minimum display			
Gram	250	0.0001	101	0.00001			
Milligram	250000	0.1	101000	0.01			
Ounce (Avoir)	8.82	0.00001	3.56	0.000001			
Troy Ounce	8.04	0.00001	3.25	0.000001			
Metric Carat	1250	0.001	505	0.0001			
Momme	66.7	0.0001	26.9	0.00001			
Pennyweight	160.8	0.0001	64.9	0.00001			
Grain (UK)	3858	0.002	1559	0.0002			
Tael (HK general, Singapore)	6.61	0.00001	2.67	0.000001			
Tael (HK jewelry)	6.68	0.00001	2.70	0.000001			
Tael (Taiwan)	6.67	0.00001	2.69	0.000001			
Tael (China)	8.00	0.00001	3.23	0.000001			
Tola (India)	21.43	0.00001	8.66	0.000001			
Messghal	53.3	0.0001	21.5	0.00001			

	GH-202					
Unit	Standar	rd range	Precision range			
Crint	Capacity	Minimum display	Capacity	Minimum display		
Gram	220	0.0001	51	0.00001		
Milligram	220000	0.1	51000	0.01		
Ounce (Avoir)	7.76	0.00001	1.80	0.000001		
Troy Ounce	7.07	0.00001	1.64	0.000001		
Metric Carat	1100	0.001	255	0.0001		
Momme	58.7	0.0001	13.6	0.00001		
Pennyweight	141.5	0.0001	32.8	0.00001		
Grain (UK)	3395	0.002	787	0.0002		
Tael (HK general, Singapore)	5.82	0.00001	1.35	0.000001		
Tael (HK jewelry)	5.88	0.00001	1.36	0.000001		
Tael (Taiwan)	5.87	0.00001	1.36	0.000001		
Tael (China)	7.04	0.00001	1.63	0.000001		
Tola (India)	18.86	0.00001	4.37	0.000001		
Messghal	46.9	0.0001	10.9	0.00001		

	GH-300	GH-200	GH-120			
Unit		Capacity				
Gram	320	220	120	0.0001		
Milligram	320000	220000	120000	0.1		
Ounce (Avoir)	11.29	7.76	4.23	0.00001		
Troy Ounce	10.29	7.07	3.86	0.00001		
Metric Carat	1600	1100	600	0.001		
Momme	85.3	58.7	32.0	0.0001		
Pennyweight	205.8	141.5	77.2	0.0001		
Grain (UK)	4938	3395	1852	0.002		
Tael (HK general, Singapore)	8.47	5.82	3.17	0.00001		
Tael (HK jewelry)	8.55	5.88	3.21	0.00001		
Tael (Taiwan)	8.53	5.87	3.20	0.00001		
Tael (China)	10.24	7.04	3.84	0.00001		
Tola (India)	27.44	18.86	10.29	0.00001		
Messghal	68.3	46.9	25.6	0.0001		

5.2. Changing the Units

The units or modes can be selected and stored in the function table. The sequence of displaying these can be arranged to fit the frequency of use.

The units stored are maintained in non-volatile memory, even if the AC adapter is removed.

- 1 Press and hold the RANGE key until bASFnc of the function table is displayed in the weighing mode, then release the key.
- 2 Press the RANGE key several times to display Unit.
- 3 Press the PRINT key to enter the unit selection mode.
- 4 Specify a unit or mode in the order to be displayed using the following keys.

RANGE key ······ To display the units sequentially.


RE-ZERO key ···· To specify a unit or mode.

The stabilization indicator o appears when the displayed unit or mode is specified.

Examples

Unit	Display	
Gram	g	°נוח ול g
Milligram	mg	°Un ıŁ mg
Counting mode	pcs	°Un ıŁ pcs
Percent mode	%	°ט יד %
Densitymode	d	°Un ıŁ d


- 5 Press the PRINT key to store the units or modes. The balance displays End and then displays the next menu item of the function table.
- 6 Press the CAL key to exit the function table. Then the balance returns to the weighing mode with the selected unit.

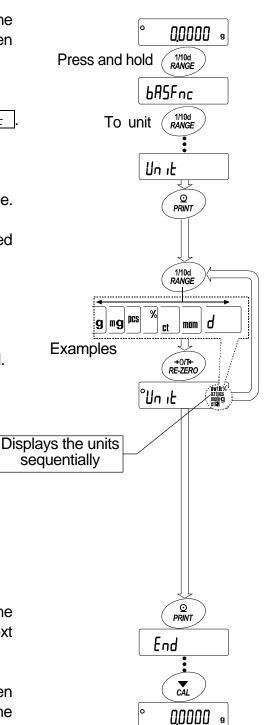

5.2.1. Unit setting example

The example below sets the units in the order with g (gram) as the first unit followed by pcs (counting mode).

- 1 Press and hold the RANGE key until bR5Fnc of the function table is displayed in the weighing mode, then release the key.
- 2 Press the RANGE key several times to display Unit
- 3 Press the PRINT key to enter the unit selection mode.
- 4 Press the RE-ZERO key to specify the unit of g.
 The stabilization indicator **O** appears when the unit is specified.
- 5 Press the RANGE key several times to display Unit PDS.
- 6 Press the RE-ZERO key to specify the unit of pcs.
 The stabilization indicator O appears when the unit is specified.
- 7 Press the PRINT key to store the units.
 The balance displays [End] and then displays the next menu item of the function table.
- 8 Press the CAL key to exit the function table. Then the balance returns to the weighing mode with g, the unit selected first.
- 9 Press the MODE key to switch between g and pcs (g→pcs).

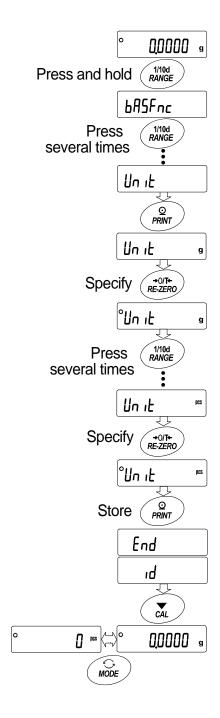

6. Weighing

Precautions for the weighing operation

- Press the RE-ZERO key each time, before placing a sample on the weighing pan, to prevent possible errors.
- Place a sample in the center of the weighing pan gently.
- Temperature changes during measurement may cause weighing error.
- Shorten the operation time as much as possible. (Opening and closing door, putting and removing sample)
- Use a pair of tweezers to avoid a temperature change due to having your hand in the weighing chamber.
- Material with an electrostatic charge or that is magnetic may cause a weighing error.
- Do not press keys with a sharp instrument (such as a pencil or ball point pen).
- Do not drop things on the pan, or place a weight on the pan that is beyond the weighing range of the balance.
- Calibrate your balance periodically to maintain weighing accuracy. Refer to section "8.
 Calibration".
- Keep the area clean and dry.
- Consider section "3. Precautions" for the weighing operation.
- For precision weighing, keep the AC adapter connected to the balance.

6.1. Basic Operation (Gram Mode)

Read section "4. Display symbols and Key operation" before operation.


Note When turning on the balance with a container placed on the pan, the tare function sets the display to zero automatically.

- 1 Turn on the balance using the ON:OFF key.
- 2 Select a preset unit (g or mg) using the MODE key.
- 3 Place the container on the weighing pan, if necessary.

Press the RE-ZERO key to cancel the weight (tare). Then zero is displayed.

Container: A vessel placed on the pan, but not to be included in the weighing data.

- 4 Place a sample on the pan or in the container.
- 5 Wait for the stabilization indicator **O** to be displayed, then read the value.
- 6 Remove the sample and container from the pan.

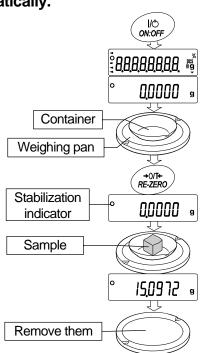


6.2. Dual Range


	Weighing range					Available minimum displa			
	GH-252		GH-202		Available	minimum	display		
Precision range	0 g	to	101 g	0 g	to	51 g	0.01 mg	0.1 mg	1 mg
Standard range	101 g	to	250 g	51 g to 220 g			0.1 mg	1 mg	

The GH-252 and GH-202 are equipped with two ranges "precision range" and "standard range".

- When weighing is started by pressing the ON:OFF key, the minimum display will be 0.1 mg.
- Pressing the RANGE key will switch these ranges alternately.

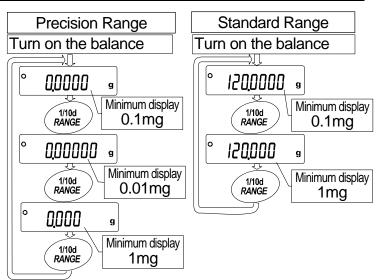

- When a sample is weighed using the precision range with a minimum display of 0.01 mg and the weight value exceeds the precision range value, the minimum display changes to 0.1 mg of the standard range. When removing the sample, the minimum display changes to 0.01 mg of the precision range automatically.
- When a tare weight (container mass value) exceeds the precision range value, even if the sample is within the precision range value, 0.01 mg of the precision range can not be selected for the minimum display. In order to select 0.01 mg of the precision range, remove the tare weight and press the RE-ZERO key to cancel it.
- When a minimum display of 0.1 mg or 1 mg is selected by the RANGE key, the minimum display is maintained even if the range is changed during weighing.

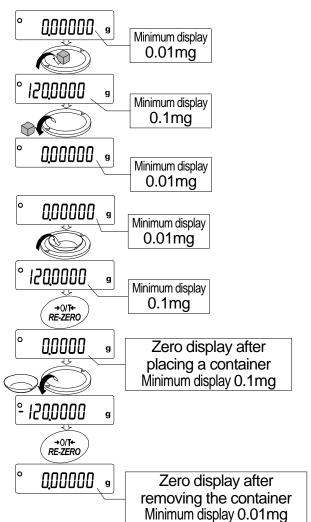

6.3. Counting Mode (PCS)

This is the mode to determine the number of objects in a sample based on the standard sample unit mass. The unit mass means an average mass of the samples. The smaller the variation in the samples, the more accurate the count will be. The balance is equipped with the Automatic Counting Accuracy Improvement (ACAI) function to improve the counting accuracy.

Notes

- Use samples with a unit mass of 1 mg or more for counting.
- If the sample unit mass variable is too large, it may cause a counting error.
- To improve the counting performance, use the ACAI function frequently or divide the samples into several groups and count each group.

Selecting the counting mode

1 Press the MODE key to select the unit pcs (counting mode).

Storing a sample unit mass (Weighing input mode)


- 2 Press the RANGE key to enter the sample unit mass storing mode.
- 3 To select the number of samples using the RANGE key. It may be set to 10, 25, 50 or 100.
 - Advise A greater number of samples will yield more accurate counting result.
- 4 Place a container on the weighing pan, if nec essary. Press the RE-ZERO key to cancel the weight (tare). The number specified in step 3 appears.
 - Example: 25 0 pcs is displayed if 25 is selected in step 3.
- 5 Place the number of samples specified on the pan. In this example, 25 pieces.
- Wait for the stabilization indicator to be displayed. Press the PRINT key to calculate and store the unit mass. Then the balance displays 25 pcs and is set to count samples with this unit mass. (The sample unit mass is stored in non-volatile memory, and is maintained even if the AC adapter is removed.) To improve the accuracy of the unit mass, proceed to step 8.

Notes

- If the balance judges that the mass of the samples is too light (under 0.0001g) and can not be stored as the unit mass, it displays La.
- If the balance judges that the mass of the samples is too light to acquire accurate weighing, it displays an error requiring the addition of more samples to the specified number.
 Example: 50 pcs appears, requiring 25 more samples. Add 25 samples and press the PRINT key. When the unit mass is stored correctly, the balance proceedes to the counting mode.

Counting operation

7 Place the samples to be counted on the pan.

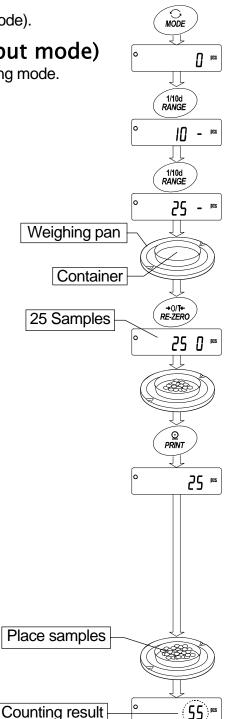


Counting mode using the ACAI function

The ACAI is a function that improves the accuracy of the unit mass automatically by increasing the number of samples as the counting process.

ACAI: Automatic Counting Accuracy Improvement

- 8 If a few more samples are added, the processing indicator turns on. To prevent an error, add three or more. The processing indicator does not turn on if overloaded. Try to add the same number of samples as displayed.
- 9 The balance re-calculates the unit mass while the processing indicator is blinking. Do not touch the balance or samples on the pan until the processing indicator turns off.
- 10 Counting accuracy is improved when the processing indicator turns off. Each time the above operation is performed, a more accurate unit mass will be obtained. There is no definite upper limit to the ACAI range for the number of samples exceeding 100. Try to add the same number of samples as displayed.
- 11 Remove all the samples used in ACAI and proceed with the counting operation using the improved unit mass.

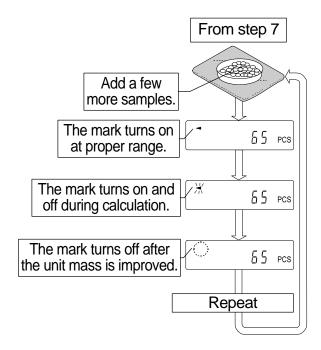

Note ACAI will not function on the unit mass entered using the keys, or digital input mode.

6.4. Percent Mode (%)

The percent mode displays the weight value in percentage compared to a 100% reference mass and is used for target weighing or checking the sample variance.

Selecting the percent mode

1 Press the MODE key to select the unit % (Percent mode). If the percent mode can not be selected, refer to "5. Weighing Units".

Storing the 100% reference mass

- 2 Press the RANGE key to enter the 100% reference mass storing mode.
 - Even in the storing mode, pressing the MODE key will switch to the next mode.
- 3 Place a container on the weighing pan, if necessary. Press the RE-ZERO key to cancel the weight (tare). The balance displays 100 0%.
- 4 Place the sample to be set as the 100% reference mass on the pan or in the container.
- 5 Press the PRINT key to store the reference mass. The balance displays 1000 %. (The decimal point position depends on the reference value. The reference mass is stored in non-volatile memory, and is maintained even if the AC adapter is removed.)

Note


Position of decimal point can be changed by 100% mass.

			<u> </u>
	100	Minimum display	
0.0100g	to	0.0999g	1%
0.1000g	to	0.9999g	0.1%
1.0000g	to	weighing capacity	0.01%

- □ If the balance judges that the mass of the sample is too light (under 0.01g) to be used as a reference, it displays □ L□ □.
- A 100% reference mass can be stored in the non-volatile memory and is maintained even if the AC adapter is removed.
- 6 Remove the sample.

Reading the percentage


7 Place a sample to be compared to the reference mass on sthe pan. The displayed percentage is based on the 100% reference mass.

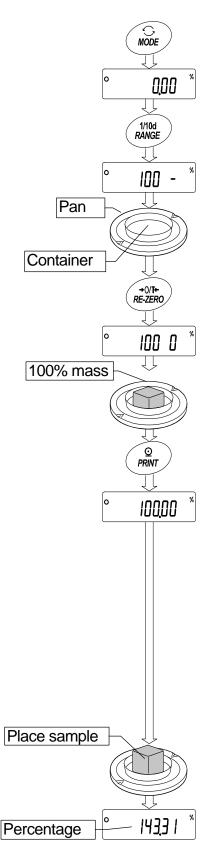

7. Response Adjustment

This function stabilizes the weight value, reducing the influence on weighing that is caused by drafts and/or vibration at the place where the balance is installed. This function adjusts by automatically analyzing the environment or by hand-operation. The function has three stages as follows:


Indicator	Parameter	Response	Stability
FAST	Cand O	Fast response,	Sensitive value
MID.	Cond I	1	•
SLOW	Cond 2	Slow response,	Stable value

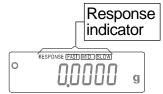

7.1. Automatic Response Adjustment

This function automatically updates the response adjustment by analyzing the influence of the environment using the internal mass.

- 1 Press and hold the MODE key until RESPONSE is displayed, then release the key.
- 2 The balance automatically sets the response characteristic.
 Caution Do not allow vibration or drafts to affect the balance during adjustment.
- 3 After automatic adjustment, the balance displays <code>End</code>, returns to the weighing mode and displays the updated response indicator for about thirty seconds.

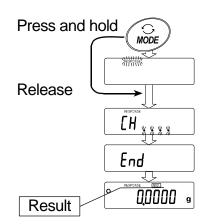

Note

- If the automatic response adjustment fails, the balance displays [[H n]]. Check the ambient conditions such as breeze and vibration, also check the weighing pan. Then, perform the adjustment again. To return to the weighing mode, press the CAL key.
- □ If there is matter on the weighing pan, the balance displays [H]. Remove them from the pan. To return to the weighing mode, press the CAL key.


Advise

If the automatic response adjustment is not helpful, try "7.2. Manual Response Adjustment".

7.2. Manual Response Adjustment

- 1 Press and hold the MODE key until RESPONSE is displayed, then release the key.

 Press the MODE key again quickly.
- 2 Select a stage of the response adjustment using the MODE key. Either FAST, MID. or SLOW can be selected.
- 3 The balance displays *End*, returns to the weighing mode and displays the updated response indicator for about thirty seconds.

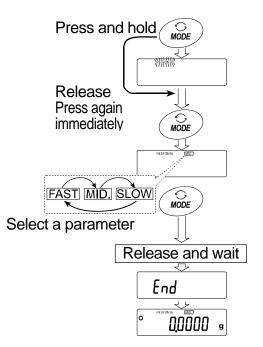

Advise

If the automatic response adjustment is not helpful, specify a parameter for "Condition (Eand) of "Environment, Display (685Fnc)" with key operation.

8. Calibration

8.1. Calibration Group

Calibration

Automatic self calibration (Calibration due to changes in temperature)

Calibration using the internal mass (One-touch calibration)

Calibration using an external weight that you have

Calibration test

Calibration test using the internal mass

Calibration test using target mass that you have

Correction of the internal mass value

Correction of the internal mass value

Caution

- Do not allow vibration or drafts to affect the balance during calibration.
- Calibration test does not perform calibration.
- □ To output the data for GLP using the RS-232C interface, set "GLP output (¬¬F□)" of "Data output (¬¬F□)". Refer to "10. Function Table". Time and date can be added to GLP report. If the time or date is not correct, adjust them. Refer to "10.7 Clock and Calendar Function".
- □ Calibration test is available only when "GLP output (יח + ם)" of "Data output (dout)" is set .
- □ The calibration and calibration test data can be stored in memory. To store them, set "Data memory (ಡਸੈ੮ਸ)". Refer to "12. Data Memory" for details.

Caution on using an external weight

- The accuracy of an external weight can influence the accuracy of weighing.
- Select a mass for calibration and calibration test from the following table.

Model	Usable calibration mass					Adjustable range
GH-120		50g,	100g*			
GH-200			100g,	200g*		-15.0 mg to +15.9 mg
GH-300			100g,	200g* ,	300g	
GH-202, GH-252	20g,	50g,	100g,	200g*		-15.00 mg to +15.99 mg

The calibration mass in **bold type**: factory setting

The calibration mass value can be adjusted within the range above.

Display

•	This indicator means in process of measuring calibration data.
	Do not allow vibration or drafts to affect the balance while the indicator is
	displayed.

About the internal mass

The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.

8.2. Automatic Self Calibration

Automatic self calibration due to changes in temperature

This function automatically calibrates the balance when the balance detects an ambient temperature change. If GLP output is selected in the function table, the balance outputs the calibration report or stores the data in memory. Automatic self calibration functions even if the display is turned off (standby state). Refer to "9-1. Permit Or Inhibit" for the operation.

Caution

- When using automatic self calibration, do not place something on the weighing pan.
- If something is on the weighing pan, the balance decides that it is in use and does not perform automatic self calibration.
- When weighing a light sample or installing the balance in a system, turn off automatic self calibration.

Note	When turning on the balance with nothing on the pan, if a sample heavier than 0.5
	g is placed on the pan, the balance detects the state that a sample is placed on the
	pan and does not perform the automatic self calibration.

_
٦

When the balance detects a change in ambient temperature, this indicator blinks and automatic self calibration is required. If the balance is not used for several minutes with this indicator blinking, the balance performs automatic self calibration. The environment will affect the time that the indicator blinks.

7				
EAL				
l – –	•	•	•	

The balance is measuring calibration data. Do not allow vibration or drafts to affect the balance while this indicator is displayed. After calibration, the balance returns to indicate the previous display.

Advise


The balance can be used while the indicator blinks. But, it is recommended that to maintain the accuracy, stop using the balance and confirm that there is nothing on the pan and allow the balance to perform self calibration.

8.3. One-Touch Calibration

This function calibrates the balance using the internal mass.

The only operation required is to press the CAL key.

- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press the CAL key to display [FIL in].
- 3 The balance performs calibration using the internal mass. Do not allow vibration or drafts to affect the balance.
- 4 The balance displays End after calibration. If the GLP output is set, the balance displays GLP and outputs the calibration report using the RS-232C interface or stores the data in memory. Refer to "GLP GLP output output (InFa)" and "Data memory (dRER)" of the function table.
- 5 The balance will automatically return to the weighing mode after calibration.
- 6 Confirm weighing accuracy using calibration test (EE 1/17).

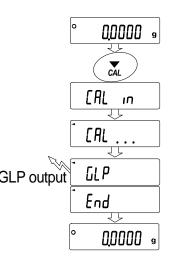


8.4. Calibration Test Using the Internal Mass

This function tests the balance accuracy using the internal mass. (Balance is not calibrated) When the GLP output is set, the calibration test report is output or stored.

- 1 Connect the AC adapter and warm up the balance at least one hour.
- 2 Press and hold the CAL key until is LE in displayed, then release the key.
- 3 The balance measures the zero point.
 Prevent vibration and drafts to affect the balance.
- 4 The measured zero point data is displayed.
- 5 The balance measures the full scale data.
 Prevent vibration and drafts to affect the balance.
- 7 The full scale data is displayed. The tolerance of the full scale data is ±0.2mg.

Model	Full scale data	
GH-120	100.000 g	
GH-200, GH-300,	200 000 a	
GH-202, GH-252	200.000 g	

- 8 When the GLP output is set, the calibration test report is output or stored.
 - □ Refer to "GLP output (¬¬F□)" and "Data memory (¬¬F□)" of the function table.
 - The zero point data and full scale data is displayed (or output) in unit of 0.0001g.
- 8 The balance automatically returns to the weighing mode.

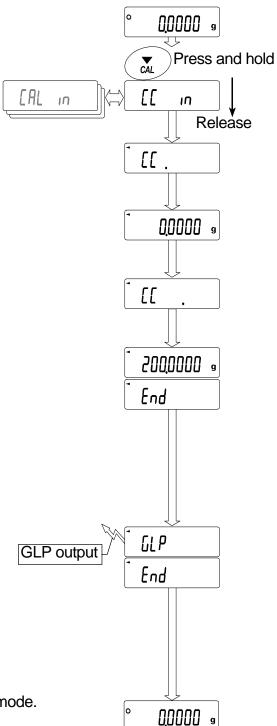

8.5. Calibration Using an External Weight

This function calibrates the balance using an external weight.

1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.

- 2 Press and hold the CAL key until [FRL out] is displayed, then release the key.
- 3 The balance displays [FRL]].
 - □ If you want to change the calibration mass, press the RANGE key and proceed to step 4.
 - If you use the calibration mass value stored in the balance, proceed to step 5.
- 4 Specify the calibration mass value as follows:

RANGE key The key to switch blinking figures.


RE-ZERO (+)key ··· The keys to select the calibration mass or MODE (-)key adjust the mass value. Refer to page 23.

PRINT key..... The key to store the new mass value. Even if the AC adapter is removed, the data is

maintained in non-volatile memory.


CAL key The key to cancel the operation and return to [FIL II].

- Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point.
 Do not allow vibration or drafts to affect the balance.
 The balance displays the calibration mass value.
- 6 Place the displayed calibration weight on the pan and press the PRINT key. The balance measures the calibration mass. Do not allow vibration or drafts to affect the balance.
- 7 The balance displays $\boxed{\mathcal{E}_{nd}}$. Remove the weight from the pan.
- 8 If the "GLP output (InFa)" the function table is set, the balance displays LLP and outputs or stores "Calibration Report". Refer to "11.2. GLP Report" for details.
- 9 The balance will automatically return to the weighing mode.
- 10 Place the calibration weight on the pan and confirm that the value displayed is correct. If it is not within the range, check the ambient conditions such as breeze and vibration also check the weighing pan. Then, repeat steps 1 to 10.

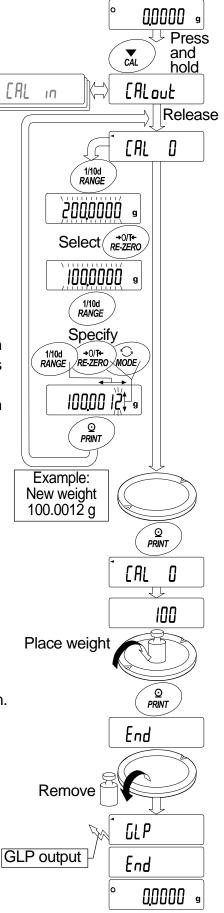

8.6. Calibration Test Using an External Weight

This function tests the weighing accuracy using an external weight. Calibration test report can be output or stored with "GLP output (IDFD)" (Calibration test does not perform calibration).

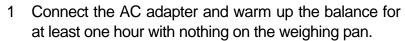

- 2 Press and hold the CAL key until [[aut is displayed, then release the key.
- 3 The balance displays [[[]].
 - □ If the target mass is changed, press the RANGE key and proceed to step 4. A list of usable weights is on page 23.
 - If current target mass value is used, proceed to step 5.
- 4 Specify the target mass value as follows:

RANGE key The key to switch blinking figures.


RE-ZERO (+)key The keys to select the target mass or adjust the mass value. Refer to page 23.

PRINT key The key to store the new mass value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.

CAL key The key to cancel the operation and return to

[[] [] []

- 5 Confirm that there is nothing on the pan and press the PRINT key. The balance measures the zero point and displays the measured value. Do not allow vibration or drafts to affect the balance. The balance displays the target mass value.
- Place the displayed target mass on the pan and press the PRINT key. The balance measures the target mass and displays the measured value. Do not allow vibration or drafts to affect the balance.
- 7 The balance displays *End*. Remove the weight from the pan.
- 8 The balance displays [GLP] and outputs or stores "calibration test report. Refer to "11.2. GLP Report" of the function table for details.
- 9 The balance will automatically return to the weighing mode.

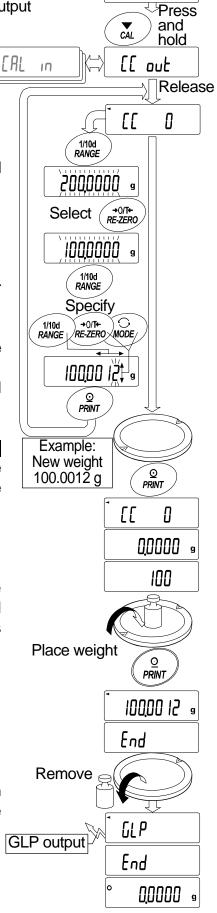

0.0000 a

8.7. Correcting the Internal Mass Value

The balance can correct the internal mass value within the range shown below. This function corrects the internal mass value to conform to an external weight. The corrected mass value is maintained in non-volatile memory even if the AC adapter is removed. The internal mass value is corrected as follows:

 Model
 Target
 Range


 GH-120
 100.000 g

 GH-200
 200.000 g

 GH-300
 200.000 g

 GH-202
 41.5 mg

 GH-252
 300.000 g

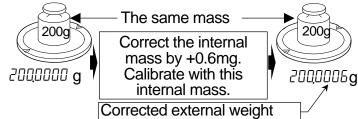

- 1 Calibrate the balance using the internal mass. (one-touch calibration)

 Example: 200.0000 g is corrected to +0.6 mg (200.0006 g). When correcting a 100 g external weight by +0.6 mg, the weight changed into 200 g, the correction value is +1.2 mg.
- 2 Press the ON:OFF key to turn off the display.
- 3 While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key. The balance displays P5.
- 4 Press the PRINT key. Then the balance displays the function switches. Set the function table switch and internal mass correction switch to " /" as shown above using the following keys.

RANGE key The key to select blinking figure.

RE-ZERO key ····· The key to change the value of the blinking figure..

PRINT key The key to store it and return to weighing mode.

CAL key The key to cancel current operation.

Function table switch

Correction switch of internal mass

5 Press and hold the RANGE key to enter the function table and release the key when basen is displayed.

- 6 Press the RANGE key several times until [5 in] is displayed, then release the key.
- 7 Press PRINT key. Correct the internal mass value using the following keys.


RE-ZERO (+)key ·· The key to select the value.

MODE (-)key ······· The key to select the value.

PRINT key The key to store the new value and display the next menu item of the function table.

CAL key.....The key to cancel this correction and display the next menu item of the function table.

- 8 Press the CAL key to return the weighing mode.
- $\,9\,\,$ Press the $\,$ CAL $\,$ key to calibrate the balance using the internal mass.
- 10 Check the correction that has been performed properly with the external weight. If the value is incorrect, repeat the correction.

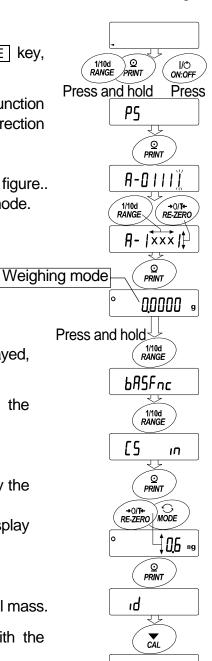

0,0000 9

9. Function Switch and Initialization

9.1. Permit or Inhibit

The balance stores parameters that must not be changed unintentionally (Example: Calibration data for accurate weighing, Data for adapting to the operating environment, Control data for the RS-232C interface). There are five switches for the purpose of protecting parameters. Each switch can select either "permit" or "inhibit". "Inhibit" protects parameters against unintentional operations.

- 1 Press the ON:OFF key to turn off the display.
- 2 While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key to display 75.
- 3 Press the PRINT key. Then the balance displays the function switches.
- 4 Specify the switches using the following keys.

RANGE keyThe key to select blinking digit.

RE-ZERO key..... The key to change the parameter for the selected switch.

To inhibit changes. (Can not be used.)

To permit changes. (Can be used.)

PRINT keyThe key to store the new parameter and return to the weighing mode.

CAL key....The key to cancel current operation and return to the weighing mode.



Function table

- \square To inhibit changes to the function table.
 - To permit changes to the function table.

Calibration using the internal mass (One-touch calibration)

- To inhibit calibration using the internal mass.
- To permit calibration using the internal mass.

Calibration using the external weight

- \square To inhibit calibration using the external weight.
- To permit calibration using the external weight.

Automatic self calibration (due to changes of ambient temperature)

- To inhibit automatic self calibration.
- To permit automatic self calibration.

Internal mass correction

- To inhibit correction.
 - To permit correction.


9.2. Initializing the Balance

This function returns the following parameters to factory settings.

- Calibration data
- Function table
- The sample unit mass value (counting mode),
 100% reference mass value (percent mode)
- □ The data that is stored in the balance using the data memory function
- External calibration weight and target mass value
- □ Function switch settings ("9.1. Permit Or Inhibit")
- Liquid density and temperature in the density mode

Note Be sure to calibrate the balance after initialization.

- 1 Press the ON:OFF key to turn off the display.
- While pressing and holding the PRINT key and the RANGE key, press the ON:OFF key to display P5.
- 3 Press the RANGE key to display [[Lr].
- 4 Press the PRINT key.
 To cancel this operation, press the CAL key.
- 5 Press the RE-ZERO key to display [[Lr [io]].
- 6 Press the PRINT key to initialize the balance.
 The balance will automatically return to the weighing mode.

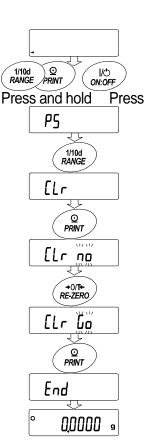

10. Function Table

The function table reads or rewrites the parameters that are stored in the balance. These parameters are maintained in non-volatile memory, even if the AC adapter is removed. The function table menu consists of two layers. The first layer is the "Class" and the second layer is the "Item".

10.1. Setting the Function Table

Display symbol and keys


0	The symbol "O" shows effective parameter.		
1/10d RANGE	When pressing and holding the key in the weighing mode, the mode enters the function table mode. The key to select the class or item in the function table mode.		
→0/T+ RE-ZERO	The key to change the parameter.		
Q PRINT	When displaying a class, the key enters an item in the class. When displaying an item, the key stores the new parameter and displays the next class.		
CAL	When displaying an item, the key cancels the new parameter and displays the next class. When displaying a class, the key exits the function table mode and returns to the weighing mode.		

Setting procedure

1	Press and hold the RANGE key until BRSFnc of the function table is displayed in the weighing mode, then release the key
2	Press the RANGE key to select a class.
3	Press the PRINT key to enter the class
4	Press the RANGE key to select a item.
5	Press the RE-ZERO key to select a parameter for the selected item.
6	If storing parameters of the selected class, press the PRINT key. Then the next class is displayed. If canceling the current operation, press the CAL key. Then the next class is displayed.
7	When specifying parameters for another class, proceed to step 2. When finishing the setting, press the CAL key to return to weighing mode.

Setting example

This example sets "Stores weighing data" for "Data memory" and "1 minute" for "Interval time".

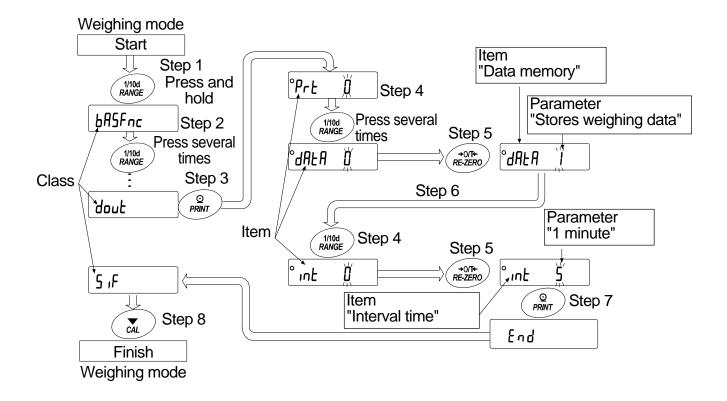

10.2. Details of the Function Table

Class	Item and Parame	ter	Desc	ription
გშ5೯იჺ Environment Display	Eand Condition	<u> </u>	Fast response, sensitive value FAST MID. Slow response, stable value SLOW	Can be changed by response adjustment.
	Stability band width	<u> </u>	Stable range is ±1 digit #1 Stable range is ±3 digits	The stabilization indicator illuminates when the display fluctuation is within the range.
	לרנ Zero tracking		OFF Normal Strong Very strong	Keeps zero display by tracking zero drift.
	5Pd Display refresh rate	- 0	5 times/second 10 times/second	Period to refresh the display
	Pnt Decimal point	- ()	Point (.) Comma (,)	Decimal point format
	P-an Auto display-ON	- 0	OFF ON	Turns on the weighing mode display when the AC adapter is connected.
EL RdJ Clock	Refer to "10.7. Clock and Calendar Function"			The time and date are added to the output data.
	Pr Ł Data output mode	- ()	Key mode	Accepts the PRINT key only when the display is stable.
		 !	Auto print mode A (Reference = zero) Auto print mode B (Reference = last stable value)	Outputs data when the display is stable and conditions of #P-P, #P-b and the reference value are met.
		3	Stream mode / Interval memory mode	With dALR D, outputs data continuously; with dALR Z, uses interval memory.
ժոսէ Data output	RP-P	• []	Plus only	Displayed value>Reference
2000	Auto print polarity	<u>-</u>	Minus only Both	Displayed value <reference displayed="" of="" regardless="" td="" value<=""></reference>
	ЯР-Ь Auto print difference	<u> </u>	10 digits 100 digits 1000 digits	Difference between reference value and displayed value #2
	ਰਸ਼ਮ Data memory	• 0 ! ?	Not used Stores weighing data Stores calibration data	Related items: Prt, int, d-no, 5-td, inFo

^{• :} Factory settings.

Example: If 1 mg display is selected using the RANGE key for the GH-300, 1 mg is one digit.

#2 Usable minimum display of the balance is one digit.

Example: In gram display, one digit is 0.01 mg for the GH-252 and 0.1 mg for the GH-300.

^{#1} The unit of minimum display is digit.

Class	Item and Parame	ter	Desc	ription
		- []	Every measurement	
			2 seconds	
		2	5 seconds	
	1	3	10 seconds	
	int	4	30 seconds	Interval time in the interval memory mode when using Prt 3, dRtR /
	Interval time	5	1 minute	Thous when doing ** E 3, Bile!
		Б	2 minute	
		7	5 minute	
		8	10 minute	
	d-no	- []	No output	Defeate #40 Dete Messes #
	Data number output		Output	Refer to "12. Data Memory"
		. []	No output	Selects whether or not the time
dout	5-Ed		Time only	
Data output	Time/Date output	2	Date only	or date is added to the weighting data. Refer to "10.7. Clock and
	•	3	Time and date	Calendar Function" for details.
	5- ıd	- []	No output	Selects whether or not the ID
	ID number output		Output	number is output.
	PUSE	- []	No pause	
	Data output pause		Pause (1.6 seconds)	Selects the data output interval.
	RL-F	- []	Not used	Selects whether or not automatic
	Auto feed		Used	feed is performed.
	ınFa	- []	No output	Selects GLP output method.
	GLP output	- 1	AD-8121 format	Refer to "11. ID Number And GLP
	GLP output	5	General data format	Report " for details.
	Ar-d	- []	Not used	Adjusts zero automatically after data
	Zero after output	1	Used	is output
		0	600 bps	
		1	1200 bps	
	6P5	• 2	2400 bps	
	Baud rate	3	4800 bps	
		4	9600 bps	
		5	19200 bps	
	66Pr	- []	7 bits, even	
5 ıF	Data bit, parity bit	1	7 bits, odd	
Serial		2	8 bits, none	
interface	[rLF	- []	CR LF	CR: ASCII code 0Dh
	Terminator	1	CR	LF: ASCII code 0Ah
		- []	A&D standard format	
			DP format	
	Ł YPE	2	KF format	Refer to "10.5. Description of Item
	LJIL			
	Data format	3	MT format	"Data Format".

■ : Factory settings.

The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number. Caution

	E-UP	0	No limit	Selects the wait time to receive a
5 ₁ F	Timeout	- /	1 second	command.
Serial	Er[d	- []	No output	AK: ASCII code 06h
interface	AK, Error code		Output	AK. ASCII CODE OOII
lilleriace	CŁ5	- []	Not used	Controls CTS and RTS.
	CTS, RTS control		Used	Controls C13 and K13.
d5 Fnc	Ldin	- []	Water temperature	Available only when density mode is
Density function	Liquid density input		Liquid density	selected
Un ıE			Refer to "5. Weighing Units".	
Unit			Neier to 3. Weighing Offics .	
[5 in				mass value correction switch is set
Internal mass of	orrection		to 1. Refer to 8. Calibration.	
ıd			Refer to "11. ID Number And GLF	D Papart"
ID number se	tting		Neiei to TT. ID Number And GEF	Nepoll.

^{• :} Factory settings. Digit is a unit of minimum weighing value.

Caution

The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

10.3. Description of the Class "Environment, Display"

Condition ([ond)

Cand () This param

1

This parameter is for sensitive response to the fluctuation of a mass value. Used for powder target mass, weighing a very light sample or when quick response weighing is required. After setting, the balance displays FAST.

Cand 2


This parameter is for stable weighing with slow response. Used to prevent a mass value from drifting due to vibration or drafts. After setting, the balance displays SLOW.

Notes In automatic response adjustment, this parameter is selected automatically.

Stability band width (5b-b)

This item controls the width to regard a mass value as a stable value. When the fluctuation per second is less than the parameter, the balance displays the stabilization indicator and outputs or stores the data. The parameter influences the "Auto print mode"

5½-b 🖟 This parameter is used for sensitive response of the stabilization indicator. Used for exact weighing.

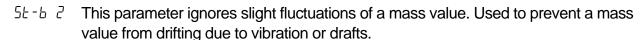

Zero tracking (*Lrc*)

This function tracks zero point drift caused by changes in the environment and stabilizes the zero point. When the weighing data is only a few digits, turn the function off for accurate weighing.

- The tracking function is not used. Used for weighing a very light sample.
- Erc / The normal tracking function is used.
- $\[\]$ The strong tracking function is used.
- *Erc* ∃ The very strong tracking function is used. Used for stable zero display.

Display refresh rate (5Pd)

The period to refresh the display. This parameter influences "Baud rate", "Data output pause" and the data output rate of "Stream mode".

Decimal point (Pnt)

The decimal point format can be selected.

Auto display-ON (P-on)

When the AC adapter is plugged in, the display is automatically turned on without the ON:OFF key operation, to display the weighing mode. Used when the balance is built into an automated system. one hour warm up is necessary for accurate weighing.

10.4. Description of the Item "Data Output Mode"

The parameter setting of "Data output mode (PrE)" applies to the performance when the "Data memory (dRER)" parameter is set to "Z" (to store the weighing data) and when the data is transmitted using the RS-232C interface.

Key mode

When the PRINT key is pressed with the stabilization indictor turned on, the balance outputs or stores the weighing data and the display blinks one time.

Required setting doub Prt (1) Key mode

Auto print modes A and B

When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and reference value are met, the balance outputs or stores the weighing data. When the PRINT key is pressed with the stabilization indictor turned on, the balance outputs or stores the data and the display blinks one time.

Auto print modes A

Example For weighing each time a sample is placed and removed, with "Rr - d"

set to " /" (to adjust zero after the data is output).

Required setting doub Prb Auto print mode A (reference = zero)

dauk RP-P Auto print polarity
dauk RP-b Auto print difference
dauk Rr-d | Zero after output

Auto print modes B

Example For weighing while a sample is added.

Required setting doub Prb 2 Auto print mode B (reference = last stable value)

dout RP-P Auto print polarity
dout RP-b Auto print difference

Stream mode

The balance outputs the weighing data continuously regardless of the display condition. When the display refresh rate is set to 5 times / second (5Pd \Box), the data output rate is also set to the same 5 times / second. The display does not blink in this mode. The interval memory mode is used when the "Data memory (dRER)" parameter is set to "!" (to store the weighing data).

Example For monitoring data on a computer.

Required setting doub Prt 3 Stream mode

לסטב dRER וו Data memory function is not used

bRSFnc SPd Display refresh rate

5 if bP5 Baud rate

Caution The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

Interval memory mode

The weighing data is periodically stored in memory.

Example For periodical weighing without a personal computer command and

to output all of the data, to a computer, at one time.

The GH series can use time and date with "Time/Date output (5-6)".

Required setting dout Prt 3 Interval memory mode

dout dRtR 2 Data memory function is used

dout וחל Interval time

Optional setting doub 5-bd/, 2, or 3 Adds the time and date.

10.5. Description of the Item "Data Format"

A&D standard format

5 if LYPE D


This format is used when the peripheral equipment can receive the A&D format.

If an AD-8121B is used, set the printer to MODE 1 or 2.

□ This format consists of fifteen or sixteen characters excluding the terminator.

When numerical characters without decimal point are exceeded eight characters for the GH-252, the format becomes sixteen characters.

- A header of two characters indicates the balance condition.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is applied.
- The unit, consisting of three characters, follows the data.

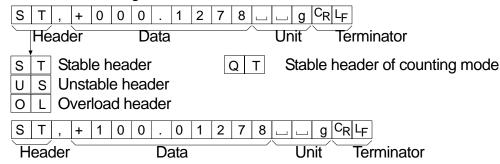


DP (Dump print) format

5 if LYPE I

This format is used when the peripheral equipment can not receive the A&D format. If an AD-8121B is used, set the printer to MODE 3.


- This format consists of sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition. No overload header is used.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- The unit, consisting of three characters, follows the data.

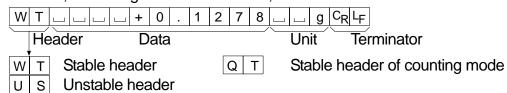

KF format 5 iF EYPE 2


This is the Karl-Fischer moisture meter format and is used when the peripheral equipment can only communicate using this format.

- This format consists of fourteen characters excluding the terminator.
- This format has no header characters.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- This format outputs the unit only for a stable value.

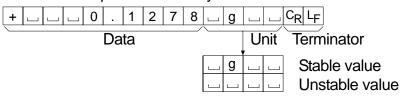

MT format 5 if EYPE 3

- A header of two characters indicates the balance condition.
- □ The polarity sign is used only for negative data.
- The weighing data uses spaces in place of the leading zeros.
- The character length of this format changes dependent upon the unit

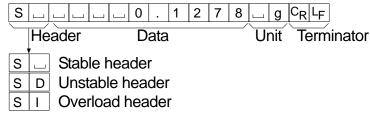

NU (numerical) format

SIF HYPE Y

This format outputs only numerical data.

- This format consists of ten characters excluding the terminator.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is used.



CSV format

SIF LYPE S

- □ This format separates the data of A&D standard format and the unit by a comma (,).
- □ This format outputs the unit even when the data is overloaded.
- □ When a comma (,) is selected for the decimal point, the separators are set to semicolon (;).
- □ When the ID number, data number, time and date are added at "Data output (dout)" of the function table, outputs ID number, data number, date, time and weighing data in this order and separates each item by a comma and treats all the items as one group of data.

LAB-0123, No,012, 2004/07/01, 12:34:56, ST,+0000.1278, g<CR><LF>
ID number Data number Date Time Weighing data

S	Т	,	+	0	0	0		1	2	7	8	,			g	C_{R}	LF	
0	L	,	+	9	9	9	9	9	9	9	Е	+	1	9	,	ш	ш	g C _R L _F

ID number

dout 5-id 1

The number to identify a specific balance.

This format consists of eight characters excluding the terminator.

Data number

dout d-no 1

This format outputs the data number just before the data is transmitted using the RS-232C interface.

- □ This format consists of six characters excluding the terminator.
- \Box When CSV format (5 ,F & BPE 5) is selected, the period (.) is replaced with a comma (,).

N o . 0 0 1 C_RL_F

Data number Terminator

Date

dout 5-td 2 or 3

□ The date output order can be changed in "Time/Date output (5-₺d)" and "Clock (₺₺ ₭₺d)". The year is output in a four-digit format.

2 0 0 4 / 0 7 / 0 1 C_RL_F

Time

dout 5-td 1 or 3

 $\hfill\Box$ This format outputs time in 24-hour format.

1 2 : 3 4 : 5 6 C_R L_F

10.6. Data Format Examples

Stable

0	0, 1278	_
	ט, וכ ום	g

A&D	S	Т	,	+	0	0	0		1	2	7	8			g	C_{R}	LF	
DP	W	Т		Ĺ			+	0		1	2	7	8]		g	C_R	L_{F}
KF	+	J			0		1	2	7	8		g			C_R	LF		
MT	S	J		L	Ĺ	Ĺ	0		1	2	7	8		g	C_R	LF		
NU	+	0	0	0	0		1	2	7	8	C_{R}	LF						

Unstable

0	-	18,3690	g
---	---	---------	---

A&D	U	S	,	-	0	1	8		3	6	9	0	ш	ш	g	C_{R}	L_{F}	
DP	U	S			Г	-	1	8		3	6	9	0]		g	C_{R}	ĿF
KF	-]	1	8		3	6	9	0]	C_R	LF		
MT	S	D]		1	1	8		3	6	9	0]	g	C_{R}	LF		
NU	-	0	0	1	8		3	6	თ	0	C_{R}	LF						

Overload


Positive error



A&D	0	L	,	+	9	9	9	9	9	9	9	Е	+	1	9	C_{R}	LF	
DP	J]	J]		J	J	Ш	J]]]]]]	C_{R}	LF
KF			J			Ĺ	Н]]]	J]	C_{R}	LF		
MT	S	ı	+	c_R	۱F													
NU	+	9	9	9	9	9	9	9	9	9	C_{R}	LБ						

Overload

Negative error

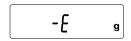

A&D DP	0	L	,	•	9	9	9	9	9	9	9	Ε	+	1	9	C_{R}	LF	
DP	ш						ш	-	Е	ш	Г	Г	ш]			C_R	LF
KF]				Г	L]		Г		ш]	C_{R}	LF		
MT	S	Ι	-	C_R	L_{F}													
NU	-	9	9	9	9	9	9	9	9	9	C_{R}	ЬF						

□ Space, ASCII 20h

^CR Carriage Return, ASCII 0Dh

Line Feed, ASCII 0Ah

Units

		A&D	D.P.	KF	MT
g	g	니니 g	니니 g	$\Box g \Box \Box$	∟ g
mg	mg	∟ m g	∟ m g	\square m g \square	_ m g
Counting mode	pcs	□ P C	⊔ P C	∟ p c s	⊔ P C S
Precent mode	%	니니%	山山%	니%니니	山%
Ounce (Avoir)	OZ	<u></u> □ 0 Z	⊔ O Z	니 O Z 니	⊔ O Z
Troy Ounce	ozt	o z t	o z t	니 o z t	니 o z t
Metric Carat	ct	_ c t	∟ c t	∟ c t ∟	∟ c t
Momme	mom	m o m	m o m	∟ m o m	∟ m o
Pennyweight	dwt	d w t	d w t	ഥd w t	ᆸᅦw t
Grain	GN	□GN	□GN	□ g r □	□GN
Tael (HK general, Singapore)	tl	山 t l	山 t l	⊔ t I s	⊔ t l
Tael (HK, jewelry)	tI	山 t l	∟ t l	山 t l h	山 t I
Tael (Taiwan)	tI	山 t l	山 t l	山 t l t	山 t I
Tael (China)	tl	∟ t I	⊔ t I	⊔ t I c	_ t I
Tola (India)	t	uu t		⊔ t O I	山 t
Messghal	m	m e s	m e s	⊔MS⊔	⊔ m
Density		□ D S	⊔ D S	_ DS_	□DS

[□] Space, ASCII 20h

10.7. Clock and Calendar Function

The balance is equipped with a clock and calendar function. When the "GLP output $(_{ID}F_{a})$ " parameter is set to " I" or "I" and the "Time/Date output $(5-E_{d})$ " parameter is set to " I", "I" or "I", the time and date are added to the output data. Set or confirm the time and date as follows:

Operation

- 1 Press and hold the RANGE key until bR5Fnc of the function table is displayed in the weighing mode, then release the key.
- 2 Press the RANGE key several times to display [[L RdJ]].
- 3 Press the PRINT key.

 The balance enters the mode to confirm or set the time and date.

Confirming the time

- 4 The current time is displayed with all the digits blinking.
 - □ When the time is correct and the date does not need to be confirmed, press the CAL key and proceed to step 8.
 - When the time is correct and the date is to be confirmed, press the
 RANGE key and proceed to step 6.
 - □ When the time is not correct and is to be changed, press the RE-ZERO key and proceed to step 5.

Setting the time

5 Set the time in 24-hour format using the following keys.

RANGE key The key to select the digits to change the value.

The selected digits blink.

RE-ZERO (-)key···· The key to increase the value by one.

MODE (+)key ······ The key to decrease the value by one.

PRINT key The key to store the new setting, display


End and proceed to step 6.

CAL key..... The key to cancel the new setting and proceed to step 6.

Confirming the date


- 6 The current date is displayed with all the digits blinking.
 - □ To change the display order of year ($\frac{1}{2}$), month ($\frac{1}{2}$) and day ($\frac{1}{2}$), press the MODE key. The date is output in the order as specified.
 - When the date is correct and the operation is to be finished, press the CAL key and proceed to step 8.
 - When the time is to be confirmed again, press the RANGE key and proceed back to step 4.
 - □ When the date is not correct and is to be changed, press the RE-ZERO key and proceed to step 7.

Note The year is expressed using a two-digit format. For example: The year 2004 is expressed as "04".

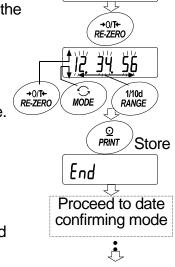

Confirm the date or

12 34 56

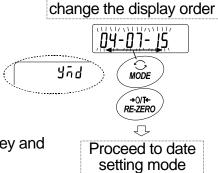

Setting the date

7 Set the date using the following keys.

RANGE key The key to select the digits to change the value.

The selected digits blink.

RE-ZERO key The key to increase the value by one.

MODE key The key to decrease the value by one.

PRINT key The key to store the new setting, display

and proceed to step 8.

CAL key The key to cancel the new setting and

proceed to step 8.

Set the date

Set the date

104-07-15

PRINT Store

End

Quitting the operation

8 The balance displays the next menu item of the function table. Press the CAL key to exit the clock and calendar function and return to the weighing mode.

Note Do not enter invalid values such as a non-existing date when setting the time and date.

When the clock backup battery has been depleted, the balance displays rtc PF. Under this condition, press any key and set the time and date. The dead battery only affects the clock and calendar function. Even so, the function works normally as long as the AC adapter is connected to the balance.


11. ID Number and GLP Report

- The ID number is used to identify the balance when Good Laboratory Practice (GLP) is used.
- □ The ID number is maintained in non-volatile memory even if the AC adapter is removed.
- The GLP output format is selected at "GLP output ($_{ID}F_{ID}$)" of the function table and can be output to a personal computer or printer using the RS-232C serial interface.
- The GLP output format includes the balance manufacturer, model, serial number, ID number, date, time and space for signature for weighing data, the weight used and results for calibration or calibration test data.
- The balance can output the following reports for GLP.
 - "Calibration report" of the calibration, using the internal mass (Calibration due to changes in temperature and one-touch calibration.)
 - "Calibration report" of the calibration, using an external weight.
 - "Calibration test report" of the calibration test, using an external weight.
 - "Title block" and "End block" for the weighing data.
- Calibration and calibration test data can be stored in memory to output several reports at the same time. Refer to "12. Data Memory" for details.
- For details on confirming and setting the time and date for the GH series. Refer to "10.7.
 Clock and Calendar Function".

11.1. Setting the ID Number



2	Press the	RANGE	key several times to display	ıd	1
_	1 1633 1116	KANGE	key several limes to display	Ü	ŀ

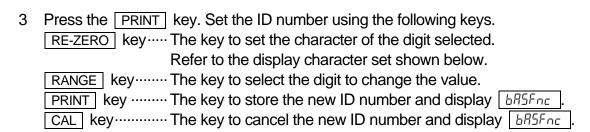

4 With b85Fnc displayed, press the CAL key to return to the weighing mode.

0	1	2	3	4	5	6	7	8	9	-	 Α	В	С	D	Ε	F	G	Н	I	J	K	L	М	N	0	Р	Q	R	S	Т	U	٧	W	Х	Υ	z
<u>[]</u>	1	2	3	4	5	5	7	8	9	-	Ŗ	Ь	Ī	덕	Ē	F	[]	Н	ı	Πļ	ዞ	Ĺ	ñ	[]	o	P	9	r	<u>י</u> ן	Ŀ	Ц	ū	<u>ا</u> ا	11	닠	Ľ

_ Space

11.2. GLP Report


Set the following parameters to output the report.

- □ Refer to "15.2. Connection to peripheral equipment" for connection to an AD-8121B.
- □ To print the report, set the "GLP output (\(\in F_B \) \) parameter to " \(\in \) and use MODE 3 of the AD-8121B. If MODE1 is used, select temporary dump print mode by pressing the \(\subseteq \text{STAT.} \) key of the AD-8121B.
- □ To output the report to a personal computer using the RS-232C interface, set the "GLP output ()" parameter to "2".
- □ If the time and date are not correct set the correct time and date in "Clock (EL Rdd)" of the function table.

Calibration report using the internal mass

Key operation

- 1 Press the CAL key to display [FRL in] and calibrate the balance automatically.
- 2 If GLP output is used, [LP] is displayed and the calibration report is output.
- 3 The balance returns to weighing mode after this calibration.

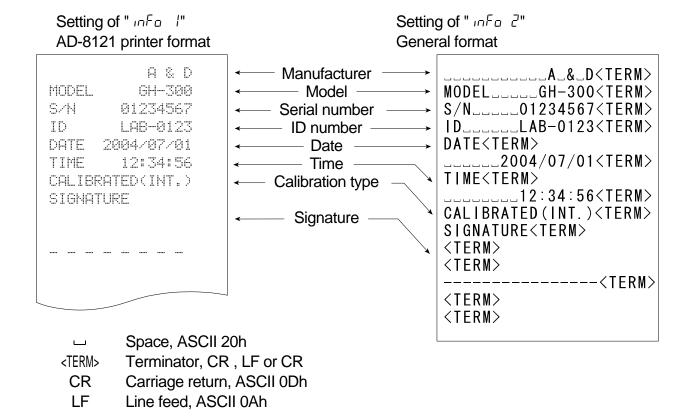

Calibration test report using an internal mass


Note Calibration test does not perform calibration.

Key operation

- 1 Press and hold the CAL key to display [[I in] and release the key.
- 2 [[] is displayed and the balance is tested automatically.
- 3 The zero point is measured and the weight value is displayed for a few seconds.
- 4 Internal mass is weighed and the weight value is displayed for a few seconds.
- 5 If GLP output is used, [LP] is displayed and the calibration test report is output.
- 6 The balance returns to weighing mode after this test.

Commad

This calibration test report can be performed with commad TST.

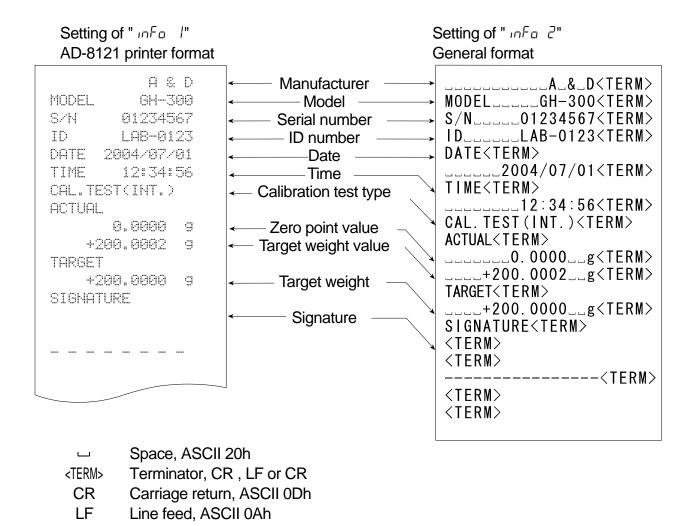

Calibration report using an external weight

Key operation

- 1 Press and hold the CAL key to display [FRL out] and release the key.
- 2 [FIL [] is displayed.
- When updating the calibration mass value, press the RANGE key and proceed to step 4. When using preset calibration mass value, proceed to step 5.
- 4 Specify calibration mass value using the following keys.


RANGE key The key to select the blinking figure

RE-ZERO (+)key ·· The key to increase the value of the blinking figure.

MODE (-)key The key to decrease the value of the blinking figure.

PRINT key The key to store the new value.

- 5 When pressing the RE-ZERO key, the zero point is measured and the weight value is displayed for a few seconds.
- 6 Place the displayed mass on the pan. Press the PRINT key to measure the mass and the weight value is displayed for a few seconds.
- 7 If GLP output is used, [LLP] is displayed and the calibration report is output.
- 8 The balance returns to weighing mode after this calibration.

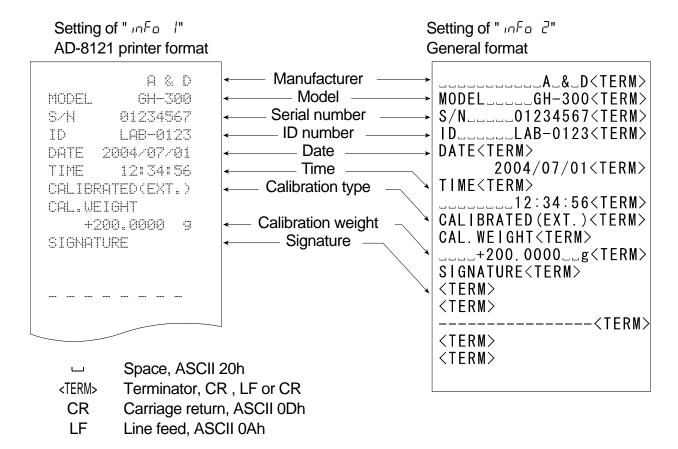

Calibration test report using an external weight

Note Calibration test does not perform calibration.

Key operation

- 1 Press and hold the CAL key to display [[out and release the key.
- 2 [FIL []] is displayed.
- When updating the target value, press the RANGE key and proceed to step 4. When using preset target value, proceed to step 5.
- 4 Specify calibration mass value using the following keys.


RANGE key The key to select the blinking figure

RE-ZERO (+)key ·· The key to increase the value of the blinking figure.

MODE (-)key The key to decrease the value of the blinking figure.

PRINT key The key to store the new value.

- 5 When pressing the RE-ZERO key, the zero point is measured and the weight value is displayed for a few seconds.
- 6 Place the displayed mass on the pan. Press the PRINT key to measure the mass and the weight value is displayed for a few seconds.
- 7 If GLP output is used, [LLP] is displayed and the calibration test report is output.
- 8 The balance returns to weighing mode after this test.

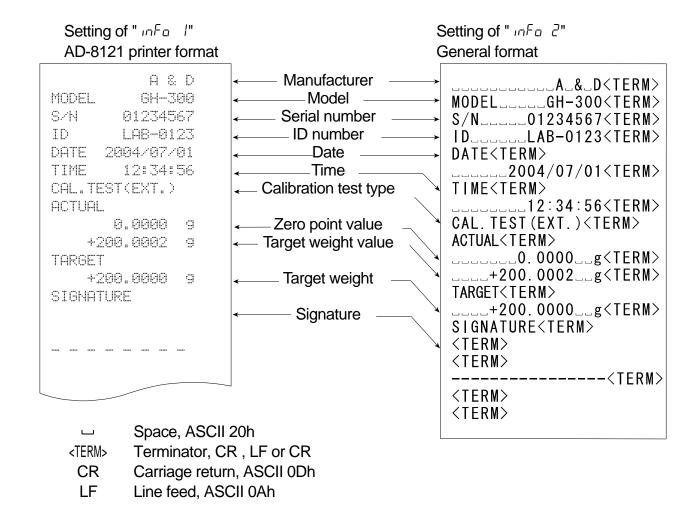

Title block and end block

When mass values are recorded as GLP data, a "Title block" is inserted at the beginning and an "End block" is inserted at the end of a group of mass values in the GLP report.


Notes

- To output the report to an AD-8121B, use MODE 3 of the AD-8121B. If MODE1 is used, select temporary dump print mode by pressing the STAT. key of the AD-8121B.
- □ If the data memory function is used (except Ⅎℛヒℛ ロ), the "Title block" and "End block" can not be output.

Key operation

- 1 With the weighing data displayed, press and hold the PRINT key, then release the key.

 5ERCE is displayed. The "Title block" is output.
- 2 The weighing data is output according to the parameter setting of the data output mode ($P_r E$) of the function table.
- 3 Press and hold the PRINT key until recent is displayed, then release the key. The "End block" is output.

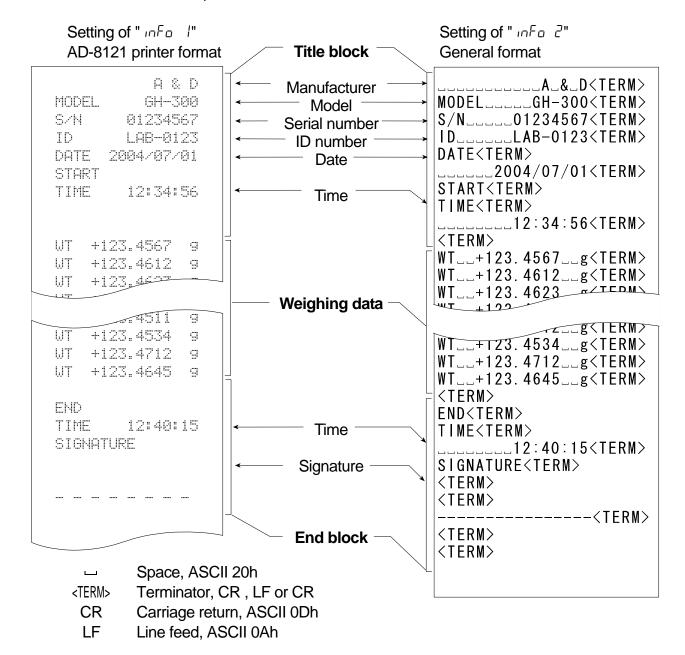

12. Data Memory

Data memory is a function to store weighing data and calibration data in memory. The data stored in memory are available for outputting at one time to a printer or personal computer.

Weighing data	Excluding date and time	Up to 200 sets
vveigi ii ig data	Including date and time	Up to 100 sets
Calibration report		
Internal calibr	ation	
External calib	ration	Last 50 sets
Calibration test rep	Lasi 50 seis	
Internal test c		
External test		

12.1. Notes on Using Data Memory

- □ To use the memory function, set the "Data memory (dRER)" parameter and "Time/Date output (5-Ed)" parameter of the function table. Refer to "10. The Function Table" for details on setting the data memory.
- □ For weighing data, the data contents to be stored and the storage capacity depend on the "Time/Date output (5-₺d)" parameter setting.

 Upper left of
- If a different type of data exists in memory, "ftr" blinks the upper left of the display. For example, you want to store weighing data but calibration data or unit mass data remains in memory.

 Under such a condition, before storing data, delete the data in memory as follows:

Releasing "[Lr" or "Err"

- 1 Press and hold the PRINT key until Ltr no with "no" blinking is displayed, then release the key.
- 2 Press the RE-ZERO key to display LLr La with "La" blinking.

 The type of data stored in memory appears in the upper left of the display as shown below:

Weighing data without time and date	-d-
Weighing data with time and date	d-F
Calibration report	H 15

- 3 Press the PRINT key to delete all the data in memory.
- 4 The balance displays $\boxed{\mathcal{E}_{nd}}$ and returns to the weighing mode.

12.2. Data Memory for Weighing Data

- The balance can store 200 sets of weighing data in memory (if time and date are added, the balance can store 100 sets). Even if the AC adapter is removed, the data is maintained in non-volatile memory.
- It is not necessary to connect the printer or personal computer to the balance continually, because the balance stores the weighing data in memory.
- The data in memory can be output at one time to a printer or personal computer.
- □ The data in memory can be displayed on the balance for confirmation.
- Data (ID number, data number, time and date) to be added to the output data can be selected in the function setting.

Storing the weighing data

Note If "[[[r"]] blinks in the upper left of the display, delete the data in memory.


- 1 Set the "Data memory (dRER)" parameter to " l".
- 2 Specify the "Time/Date output (5-৮d)" parameter whether time and date is to added or not.
- 3 The storing mode depends on the "Data output mode ($P_{\Gamma} E$)" parameter setting. Four types of storing modes are available to store data.

Key mode When the PRINT key is pressed and the displayed value is stable, the balance stores the weighing data.

Auto print modes A When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and zero point (reference value) are met, the balance stores the weighing data.

Auto print modes B When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and last stable data (reference value) are met, the balance stores the weighing data.

Interval memory mode···· Weighing data is stored at an interval specified in "Interval time (not)". Press the PRINT key to start and stop this mode.

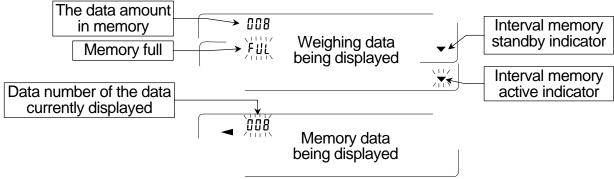

Caution

- When weighing data is being stored in memory, the data can not be output to a personal computer using the RS-232C interface.
- "FUL" indicates that memory is full or the memory capacity has been reached. More data can not be stored unless the memory data is deleted.

- Automatic self calibration can not be used while the interval memory mode is active.
- The following commands can not be used during data storage.
 - Q Query command for weighing data.
 - S Query command for stable weighing data.
 - SI Query command for weighing data.
 - SIR Query command for continuous weighing data.

Setting the function table

Parameter settings for each output mode are as follows:

Item	Data output	Auto print polarity,	Data memory	Interval time
Mode	mode	difference	function	iiileivai liiile
Key mode	Prt O	Not used	dRER I	
Auto print mode A	Prt 1	RP-R 0-2	dRER I	Not used
Auto print mode B	Prt 2	RP-6 0-2	dRER I	
Interval memory mode	Prt 3	Not used	dRER I	int 0-8

Additional parameter settings, as follows:

Data number	No	d-na O		
Data number	Yes	d-no l		
ID number	No	5- ıd 0		
number	Yes	5- ıd		

	No	5-Ed 0
Time and date	Time only	5-Ed
Time and date	Date only	5-Ed 2
	Both	5-Ed 3

Enabling the data memory function

- 1 Press and hold the RANGE key until bR5Fnc is displayed, then release the key.
- 2 Press the RANGE key several times to display doub.
- 3 Press the PRINT key.
- 4 Press the RANGE key three times to display dRLR [].
- 5 Press the RE-ZERO key to display dRER 1.
- 6 Press the PRINT key to store the setting.
- 7 Press the CAL key to return to the weighing mode.

Recalling the memory data

Confirm that the "Data memory (dRER)" parameter is set to " !".

- 1 Press and hold the PRINT key until FEERLL is displayed, then release the key.
- 2 Press the PRINT key to enter the memory recall mode. The type of data appears in the upper left of the display as shown to the right. Recall the data in memory using the following keys.

RE-ZERO key····· To proceed to the next data set.

MODE key To go back to the previous data set.

PRINT key To transmit the current data using the RS-232C interface.

With RANGE held down, press the CAL key
To delete the current data.

CAL key.....To exit the memory recall mode.

3 Press the CAL key to return to the weighing mode.

Indicators -d Weighing data without time and date

Weighing data with time and date

d-E

Transmitting all memory data at one time

Confirm that the "Serial interface (5 ,F)" parameters are set properly. Refer to "10. Function Table" and "15.2. Connection To Peripheral Equipment".

- 1 Press and hold the PRINT key until release the key.
- 2 Press the RANGE key to display aut.
- 3 Press the PRINT key to display aut no with "no" blinking.
- 5 Press the RE-ZERO key to display συξ τως with "τω" blinking.
- 6 Press the PRINT key to transmit all data using the RS-232C interface.
- 7 The balance displays [LERr] when all data is transmitted. Press the CAL key to return to the weighing mode.

Indicators -d Weighing data without time and date

Weighing data with time and date

Deleting all memory data at one time

- 1 Press and hold the PRINT key until FEERLL is displayed, then release the key.
- 2 Press the RANGE key several times to display [[LERr].
- 3 Press the PRINT key to display [[Lr ng]] with "na" blinking.
- 4 Press the RE-ZERO key to display [[L- []] with "[a" blinking.
- 5 Press the PRINT key to delete all data
- 6 The balance displays $\boxed{\textit{End}}$ and returns to the weighing mode.

12.3. Data Memory for Calibration and Calibration Test

- Calibration data (when and how it is performed) and calibration test data can be stored in memory.
- All the data in memory is available to be output at one time to a printer or personal computer.
- Up to 50 data sets of the latest calibration or calibration test can be stored. When the memory capacity has been reached, "FUL" illuminates in the upper left of the display as shown below.

Indicator FUL

Storing the calibration and calibration test data

Note If "[Lr" appears blinking in the upper left of the display, delete the data in memory.

Store the calibration and calibration test data as follows:

- 1 Set the "Data memory (♂RER)" parameter to "?".
- 2 Set the "GLP output (¬¬F□)" parameter to " /" or "?".
- With the settings above, each time calibration or calibration test is performed, the data is stored automatically.

Transmitting the memory data Note

- □ Confirm that the "Serial interface (5 ,F)" parameters are set properly.
 Refer to "10. Function Table" and "15.2. Connection To Peripheral Equipment".
- □ Confirm that the "Data memory (dREA)" parameter is set to "2".
- 1 Press and hold the PRINT key until aut is displayed, then release the key.

In the upper left of the display

H 15

- 2 Press the PRINT key to display [out no] with "no" blinking.
- 3 Press the RE-ZERO key to display aut up with "aa" blinking.
- 4 Press the PRINT key to transmit all memory data using the RS-232C interface.
- 5 The balance displays [LEAr] when all memory data is output. Press the [CAL] key to return to the weighing mode.

Deleting data stored in memory

- 1 Press and hold the PRINT key until out is displayed, then release the key.
- 2 Press the SAMPLE key to display [[LERr].

In the upper left of the display

3 Press the PRINT key to display [Lr ng with "no" blinking.

H 15


- 4 Press the RE-ZERO key to display [[Lr [[a]]]] with "[a]" blinking.
- 5 Press the PRINT key to delete all data.
- 6 The balance displays and returns to the weighing mode when all the data has been deleted.

13. Underhook


The underhook can be used for magnetic materials or density measurement. The built-in underhook is revealed by removing the cap on the bottom of the balance. Use the underhook as shown below.

Caution

- Do not apply excessive force to the underhook.
- When not in use, attach the cap to prevent dust from getting into the balance.
- Do not push the underhook upward
- When turning the balance over, the weighing pan, pan support, breeze break ring and dust plate will fall off. Remove them first.
- When the cap is removed, a weighing error may occur, because of drafts entering into the internal portion of the balance. Arrange the room condition.

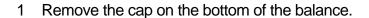

2 Hang your weighing pan on the underhook. Place the balance on the solid table.

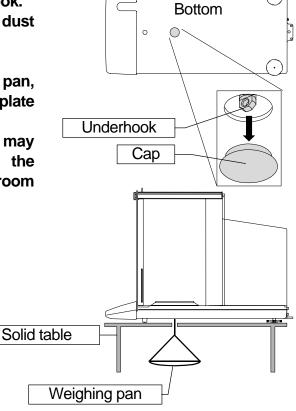

14. Density Measurement

The balance is equipped with a density mode. It calculates the density of a solid using the mass value of a sample in air and the mass value in liquid.

Note

- The density mode was not selected for use when the balance was shipped from the factory.
 To use the mode, change the function table and activate the density mode.
- □ When the density mode is selected, the response adjustment function can not be used.
- Minimum display is 0.0001 g while density mode.

Formula to obtain the density

The density can be obtained by the following formula.

 $\rho = \frac{A}{A-B} \times \rho_0$ Where ρ : Density of a sample A : Mass value of a sample in air <math display="block">B : Mass value of a sample in liquid

 ρ_0 : Density of a liquid

Prior to measurement: Changing the function table

Prior to measurement, change the function table as follows:

1 Selecting the density mode.

The density mode is available as one of the weighing units.

Select it by pressing the MODE key. To use the mode, select it (unit of Unit d.) in the function table. Refer to "5.2. Changing Units".

2 Selecting a way to set the density of a liquid. Select the liquid density input method from the function table below.

Note

The function table is available only when the density mode (d5Fnc) is selected. d5Fnc is displayed next 5 ildot F when density mode is active at Un ildot E.

Class	Item and Parameter		Description
d5 Fnc	Ldin	- []	Water temperature
Density function	Liquid density input	1	Liquid density

• is factory setting.

Entering the density of a liquid

Two ways to set the density of a liquid are available in the function table, "Liquid density input (Ld n)): by entering the water temperature or by entering the density directly.

1 Press the MODE key as necessary to select the density mode. When the density mode begins, "g (gram)" is displayed and the processing indicator (◄) blinks.



Entering the water temperature (Ld in 1)

2 The water temperature currently set (unit: °C, factory setting : 25°C) is displayed. Use the following keys to change the value.



RE-ZERO (+)key ·· The key to increase the temperature by one degree. (0°C is displayed after 99°C)

MODE (-)key The key to decrease the temperature by one degree. (99°C is displayed after 0°C)

PRINT key The key to store new water temperature, display End and return to the density mode. Proceed to Step 1.

CAL key..... The key to cancel the change and return to the density mode. Proceed to Step 1.

The relation between the water temperature and density

°C	+0	+1	+2	+3	+4	+5	+6	+7	+8	+9
0	0.99984	0.99990	0.99994	0.99996	0.99997	0.99996	0.99994	0.99990	0.99985	0.99978
10	0.99970	0.99961	0.99949	0.99938	0.99924	0.99910	0.99894	0.99877	0.99860	0.99841
20	0.99820	0.99799	0.99777	0.99754	0.99730	0.99704	0.99678	0.99651	0.99623	0.99594
30	0.99565	0.99534	0.99503	0.99470	0.99437	0.99403	0.99368	0.99333	0.99297	0.99259
40	0.99222	0.99183	0.99144	0.99104	0.99063	0.99021	0.98979	0.98936	0.98893	0.98849
50	0.98804	0.98758	0.98712	0.98665	0.98618	0.98570	0.98521	0.98471	0.98422	0.98371
60	0.98320	0.98268	0.98216	0.98163	0.98110	0.98055	0.98001	0.97946	0.97890	0.97834
70	0.97777	0.97720	0.97662	0.97603	0.97544	0.97485	0.97425	0.97364	0.97303	0.97242
80	0.97180	0.97117	0.97054	0.96991	0.96927	0.96862	0.96797	0.96731	0.96665	0.96600
90	0.96532	0.96465	0.96397	0.96328	0.96259	0.96190	0.96120	0.96050	0.95979	0.95906

Entering the density directly (Ld in 1)

The density currently set (unit: g/cm³, factory setting: 1.0000g/cm³) is displayed.

Use the following keys to change the value.

The range to set the density is 0.0000g / cm³ to 1.9999g / cm³.

9)(0000

RE-ZERO (+)key. The key to set the value of the digit selected.

RANGE (-)key······ The key to select the digit to change the value.

PRINT key The key to store the change, display End and return to the density mode. Proceed to Step 1.

CAL key..... The key to cancel the change and return to the density mode. Proceed to Step 1.

Measuring the density

Note If the liquid temperature or the type of liquid is changed during measurement, input again the value of the liquid density as necessary.

Density display is four decimal places. Minimum display can not change with the RANGE key.

Denisty is displayed after "Mass measurement in air" and "Mass measurement in liquid".

The procedure of each measurements is as follows:

- 1 Enter the density mode that "g (gram)" is displayed and the processing indicator (◄) blinks. Place nothing on both pan and press the RE-ZERO key to display zero.
- Place the sample on the pan in air. If the weight value is stored or output, press the PRINT key to store it after a stable weight value is displayed.

Press the RANGE key to decide the weight value in air and proceed to next step.

Note If negative value or E (out of range) is displayed, the RANGE key is inactive.

3 Move the sample to the pan in liquid.

If the weight value is stored or output, press the PRINT key to store it after a stable weight value is displayed.


Press the RANGE key to decide the weight value in liquid and proceed to next step.

Note If E (out of range) is displayed, the RANGE key is inactive.

4 If the density value is stored or output, press the PRINT key to store it.

Press the RANGE key to measure other sample and proceed to step 2.

- 5 If the liquid temperature or the type of liquid is changed during measurement, input again the value of the liquid density as necessary.
- 6 Press the MODE key to proceed to other modes.

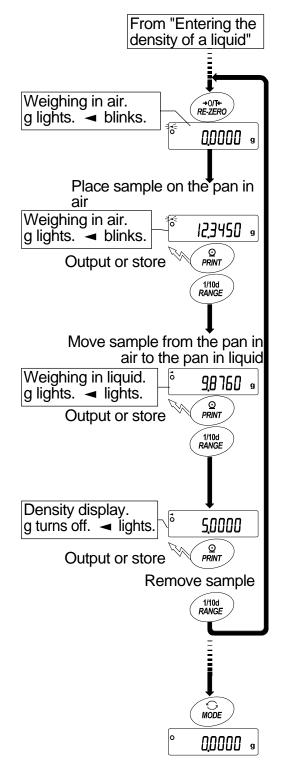

15. Standard Input and Output Interface

15.1. RS-232C Interface

The balance is a DCE device. Connect the balance to a personal computer (DTE) using a straight through cable.

Transmission system : EIA RS-232C

Transmission form : Asynchronous, bi-directional, half duplex

Transmission rate : 10 times/second or 5 times/second (same as data refresh rate)

Data format : Baud rate : 600, 1200, 2400, 4800, 9600, 19200bps

Data bits : 7 or 8 bits

Parity : Even, Odd (Data bits 7 bits)

None (Data bits 8 bits)

Stop bit : 1 bit

Code : ASCII RS-232C

1 -5V to -15V

Data bits

Start bit

RS-232C


1 -5V to -15V

Stop bit


Parity bit

D-Sub 25 pin assignments

Pin No.	Signal name	Direction	Description
1	ı	_	No connection
2	TXD	Output	Transmit data
3	RXD	Input	Receive data
4	ı	_	No connection
5	SG	-	Signal ground
6	DSR	Output	Data set ready
7	RTS	Input	Request to send
8	CTS	Output	Clear to send
9	- 1	_	No connection

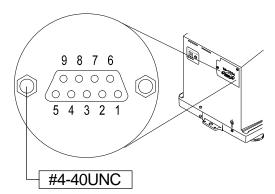

Signal names of the balance side are the same as the DTE side with TXD and RXD reversed.

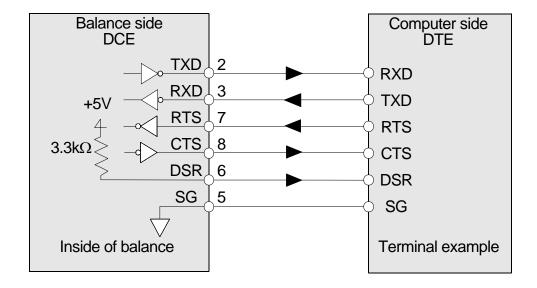

15.2. Connection to Peripheral Equipment

Connection to an AD-8121B printer

Preset the following parameters to use the AD-8121B printer.

Class	Item and Parameter	Factory settings	AD-8121B MODE 1	AD-8121B MODE 2	AD-8121B MODE 3
	Pr Ł Data output mode	0	0, 1,2	3	0, 1,2
	RP-P Auto print polarity	0	ща	Not	ш.
	RP-ь Auto print difference	1	#1	necessary	#1
dout	d-na Data number output	0	0	0	0, 1
Data output	5-Łd Time/Date output	0	0	0	0,1,2,3
	5- id ID number output	0	0	0	0, 1
	PUSE Data output pause	0	0	0	0,1 #2
	RE-F Auto feed	0	0	0	0, 1
	եր <u>Տ</u> Baud rate	2	2	2	2
5 iF	b b c Data bit, parity bit	0	0	0	0
Serial interface	ErLF Terminator	0	0	0	0
	EYPE Data format	0	0	0	1
	CTS, RTS control	0	0	0	0

^{#1} Set parameters when auto print mode A or B ($Pr \vdash l$ or 2) is selected.

Notes

□ Refer to "11.2. GLP Report" concerning print samples.

Settings of AD-8121B

50ttingo 017tb 0121b					
MODE	AD-8121B DIP switch	Description			
MODE 1	MODE \$ \$ \$	Print at receiving data. Standard mode, statistic mode			
MODE 2	MODE	Print by DATA key operation or built-in timer. Standard mode, interval mode, chart mode			
MODE 3	D D S O D O D O D O D O D O D O D O D O	Print at receiving data. Dump print mode			

DIP switch No.3: Handling unstable data

ON Print

OFF Not printed

DIP switch No.4 : Data input specifications (Interface selection)

ON Current loop

OFF RS-232C





^{#2} Set / when multiple lines are printed. Example: When appending ID number, set /.

Connection to a computer and the use of WinCT

The balance is of the DCE type (Data Communication Equipment), which can be connected to a personal computer using the RS-232C interface. Before connection, read the personal computer manual thoroughly. Use a standard DCE cable for connection (cable type: straight-through).

Using Windows Communication Tools Software (WinCT)

When Windows is used as an operating system in a personal computer, the provided WinCT software can be used to transmit the weighing data to the personal computer.

The WinCT software has two communication methods: "RsCom" and "RsKey". Refer to the WinCT instruction manual.

The current version of the WinCT can be downloaded from the A&D website.

RsCom

- RsCom can transmit commands to control the balance.
- RsCom can make bi-directional communication between the balance and a personal computer using the RS-232C interface.
- RsCom can display or store the data using a text file format. RsCom can also print the data using a printer connected to the personal computer.
- When several ports of a personal computer have balances connected, the personal computer can communicate with each balance simultaneously.
- RsCom can share a personal computer with other application software.
- RsCom can receive the balance GLP report.

RsKey

- RsKey can transmit the weighing data output from the balance directly to other application software such as Microsoft Excel.
- RsKey can be used with most application software.
- RsKey can receive the balance GLP report.

Using the WinCT software, the balance can do the following:

Analyzing the weighing data and the statistics input by "RsKey"

The weighing data can be input directly into an Excel worksheet. Then, Excel can analyze the data to obtain total, average, standard deviation, maximum and minimum value, and display them in a graph.

Controlling the balance using commands from a personal computer

By using "RsCom", the personal computer sends commands such as "re-zero" or "send weighing data" to the balance and controls the balance.

Printing the balance GLP report using your printer

The balance GLP report can be printed using a printer connected to the personal computer.

Receiving weighing data at a certain interval

The weighing data can be received at a certain interval and data characteristic with elapsed time can be obtained.

Using the balance memory function

The data can be stored in the balance's memory. Of the data stored, the weighing data and calibration data can be transmitted to a personal computer at one time.

Using a personal computer as an external indicator

With the "RsKey" test mode function, a personal computer can be used as an external weight indicator for the balance. (To do this, set the balance data output mode to stream mode.)

15.3. Commands

15.3.1. Command List

Note A command has a terminator added, that is specified using "5 if Erlf" of the function table, and is sent to the balance.

Commands to query weighing data				
С	Cancels the S or SIR command.			
Q	Requests the weighing data immediately.			
S	Requests the weighing data when stabilized.			
SI	Requests the weighing data immediately.			
SIR	Requests the weighing data continuously.			

Commands to control the balance			
CAL	Same as the CAL key.		
OFF	Turns the display off.		
ON	Turns the display on.		
P	Same as the ON:OFF key		
PRT	Same as the PRINT key		
R	Same as the RE-ZERO key (Note 1)		
RNG	Same as the RANGE key.		
TR	Tares the balance (Cancels the container's weight.) (Note 1)		
TST	Perform calibration test.		
U	Same as the MODE key		

Commands to request data		
?ID	Requests the identification number.	
?PT	Outputs the tare value. (Note 1)	
?SN	Request the serial number of the balance.	
?TN	Request the model name of the balance.	

Command to set data		
PT:***.***	Sets the tare value. (Note 1)	

 $^{\text{(Note 1)}} \; {\scriptstyle \square} \; R$ command assumes the point as zero and sets the display to zero.

- TR command cancels the tare weight when the displayed value is greater than zero.
 The canceled tare value can be read using the ?PT command.
- PT: command sets the tare value digitally.

An example of this command sets a negative target value and fills with the sample until the display becomes zero.

Step 1 Place a container.

Step 2 Set the display to zero using the R command.

Step 3 Set a target value using the PT: command.

For details, refer to page 69.

15.3.2. Acknowledge Code and Error Codes

When the "Serial interface function $(5 \ F)$ " parameter is set to " $ErEd \ I$ ", the balance outputs <AK> code or error code to each command as follows:

<AK> (06h) Acknowledge in ASCII code.

- When the balance receives a command to request data and can not process it, the balance transmits an error code (EC, Exx).
 - When the balance receives a command to request data and can process it, the balance outputs the data.
- □ When the balance receives a command to control the balance and can not process it, the balance transmits an error code (EC, Exx).
 - When the balance receives a command to control the balance and can process it, the balance transmits the acknowledge code.

Among commands to control the balance, the following transmit the acknowledge code both when the balance receives the command and when the balance has accomplished the command. If the command can not be processed properly, the balance transmits an error code (EC, Exx). This error can be released using the CAL command.

CAL command (Calibration command using internal mass)

ON command (Display ON command)

P command (Display ON/OFF command)

R command (RE-ZERO command)

TR command (Tare command)

TST command (Calibration test command)

When a communication error has occurred due to external noise, or a parity error has occurred due to transmission error, the balance transmits an error code. In this case, send the command again.

xx is error code number.

15.3.3. Control Using CTS and RTS

Depending on the " \mathcal{L} 5" parameter of "Serial interface (5 \mathcal{F})", the balance performs as follows: \mathcal{L} 5 \mathcal{D}

Regardless of whether the balance can receive a command or not, the balance keeps the CTS line HI. The balance outputs data regardless of the condition of the RTS line.

[ES 1

The CTS line is kept Hi normally. When the balance can not receive the next command (Example: while the balance is processing the last command), the balance sets the CTS line to Lo. The balance confirms the level of the RTS line before outputting a set of data. If the RTS level is Hi, the balance outputs data. If the RTS level is Lo, data is not output (The data is canceled).


15.3.4. Settings Related to RS-232C

Concerning the RS-232C, the balance has two functions: "Data output (daut)" and "Serial interface (5 ,F)". Set each function as necessary.


15.3.5. Command Examples

This example uses the "ErEd I" of "5 IF" so that the <AK> (06h) code is output.


ON command (turning on the balance)

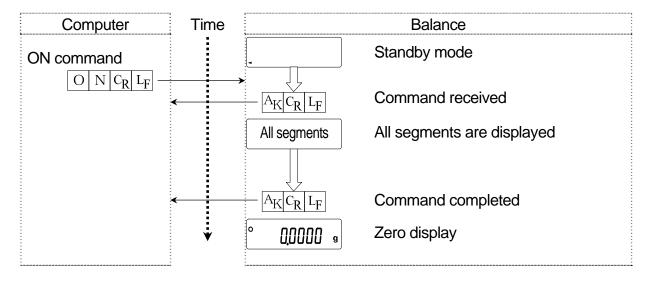

R command (Re-zeroing the display)

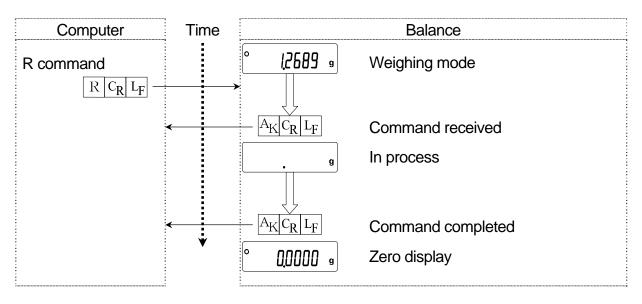


CAL command (Calibration with internal mass)

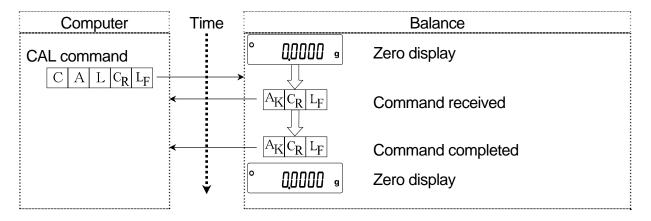


Error code

This example is of an error using the R command. "Er[d]!" is used. The balance transmits an error code when the received command can not be achieved.

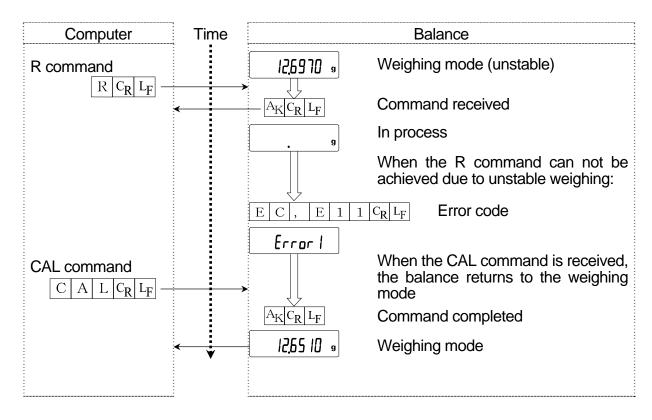


Weighing with a tare

This example uses "Er[d /" of "5 /F" so that the <AK> (06h) code is output.

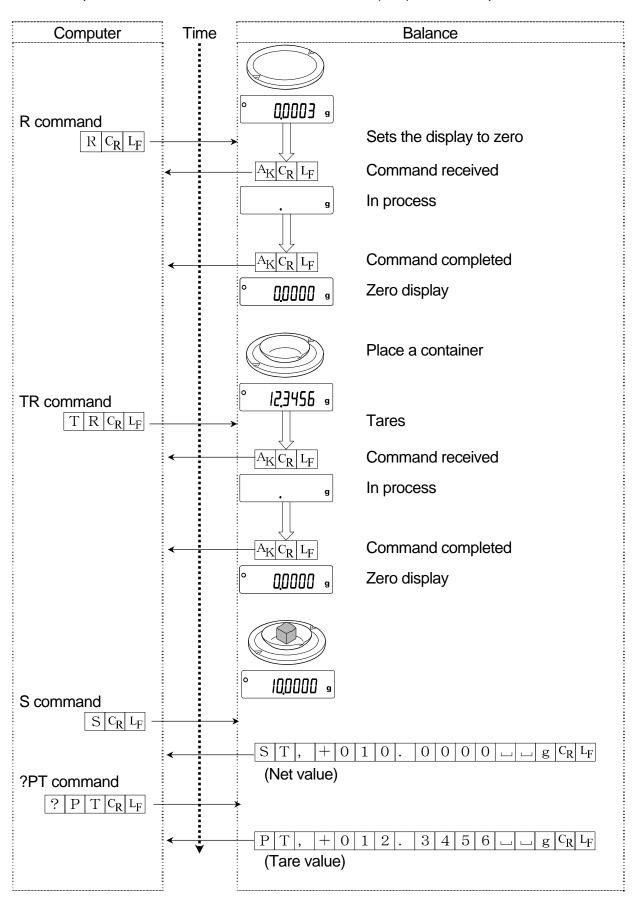

Setting a negative target value and filling with a sample until the display becomes zero

This example uses "ErEd I" of "5 F" so that the <AK> (06h) code is output.

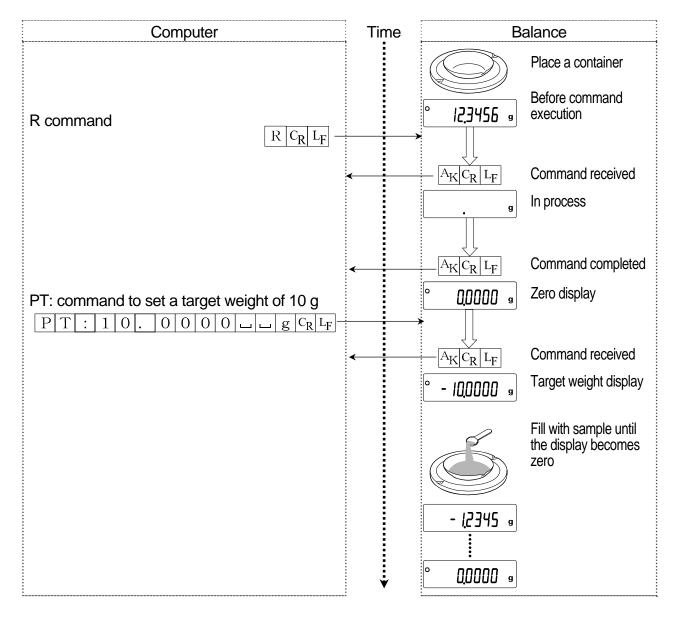

16. Maintenance

16.1. Treatment of the Balance

- Clean the balance with a lint free cloth that is moistened with warm water and a mild detergent.
- Do not use organic solvents to clean the balance.
- Do not disassemble the balance. Contact the local A&D dealer if the balance needs service or repair.
- Use the original packing material for transportation.
- Consider "3. Precautions" when operation the balance.

16.2. Error Codes

Display	Error code	Description
		Stability error
Error I	EC,E11	The balance can not stabilize due to an environmental problem. Check around the pan. Prevent vibration, drafts, temperature changes, static electricity and magnetic fields, from influencing the balance. To return to the weighing mode, press the CAL key.
		Internal mass error
Error6	EC,E16	Applying the internal mass does not yield a change in the mass value as specified. Confirm that there is nothing on the pan and perform the weighing operation from the beginning again.
		Internal mass error
[Error]	EC,E17	The internal mass application mechanism does not function properly. Perform the weighing operation from the beginning again.
		Calibration weight error
CAL E	EC,E20	The calibration weight is too heavy. Confirm the calibration mass value. Press the CAL key to return to the weighing mode.
		Calibration weight error
-CAL E	EC,E21	The calibration weight is too light. Confirm the calibration mass value. Press the CAL key to return to the weighing mode.
		Overload error
E		A sample beyond the balance weighing capacity has been placed on the pan. Remove the sample from the pan.
		Weighing pan Error
-E		The mass value is too light. Confirm that the weighing pan is properly installed and calibrate the balance.

Display Error code	Description
	Sample mass error
	The balance can not store the sample for the counting mode
Lo	or for the percent mode because it is too light.
	Use a larger sample.
	Unit mass error
	The sample unit mass for the counting mode is too light.
	Storing and using it for counting will cause a counting error.
50 - ^{pcs}	Add samples to reach the specified number and press the
188	PRINT key. Pressing the PRINT key without adding
100 - pcs	samples will shift the balance to the counting mode. But, for
	accurate counting, be sure to add samples.
	Automatic response adjustment zero error
	The automatic response adjustment can not be performed
<u> </u>	because there is something on the pan. Clear the pan. Press
	the CAL key to return to the weighing mode.
	Automatic response adjustment unstable error
<u> </u>	The automatic response adjustment can not be performed
	because the mass value is unstable.
(Check NG)	Check the ambient conditions such as breeze, vibration and
(Check 112)	magnetic fields, also check the weighing pan.
	Press the CAL key to return to the weighing mode.
	Clock battery error
	The clock backup battery has been depleted. Press any key and set the time and date. The clock and
rtc PF	calendar function works normally as long as the AC adapter
	is connected to the balance. If this error appears frequently,
	contact the local A&D dealer.
	Memory full
<u> </u>	The amount of weighing data in memory has reached the
(Blinking)	maximum capacity. Delete data in memory to store new data.
(),	For details, refer to "12. Data Memory".
	Memory full
	The amount of calibration or calibration test data in memory
FUL (Illuminated)	has reached the maximum capacity (50 sets).
(Illuminated)	The data in memory will be deleted automatically to store new
	data. For details, refer to "12. Data Memory".
	Memory type error
	Type of memory set in the function table and type of data
A. A. in	stored are different. For details, refer to "12. Data Memory".
	Memory data error
	The stored data were destroyed. To use the memory i
E.C.C.	The stored data were destroyed. To use the memory function, erase all the old data and release the error. For

Display	Error code	Description
	EC,E00	Communications error A protocol error occurred in communications. Confirm the format, baud rate and parity.
	EC,E01	Undefined command error An undefined command was received. Confirm the command.
	EC,E02	Not ready A received command can not be processed. Example: □ The balance received a "Q" command, but not in the weighing mode. □ The balance received a "Q" command while processing a RE-ZERO command. Adjust the delay time to transmit a command.
	EC,E03	Timeout error If the timeout parameter is set to "Ł-UP I", the balance did not receive the next character of a command within the time limit of one second. Confirm the communication.
	EC,E04	Excess characters error
	EC,E06	Format error A command includes incorrect data. Example: The data is numerically incorrect. Confirm the command.
	EC,E07	Parameter setting error The received data exceeds the range that the balance can accept. Confirm the parameter range of the command.
Other error code		If an error described above can not be released or other errors are displayed, contact the local A&D dealer.

16.3. Other Display



When this indicator blinks, automatic self calibration is required. The indicator blinks when the balance detects changes in ambient temperature. If the balance is not used for several minutes with this indicator blinking, the balance performs automatic self calibration. The blinking period depends on the operating environment.

Advise

The balance can be used while this indicator is blinking. We recommend that you perform automatic self calibration for precision weighing.

16.4. Checking the Balance Performance and Environment

The balance is a precision instrument. When the operating environment or the operating method is inadequate, correct weighing can not be performed. Place a sample on the pan and remove it, and repeat this several times. If the balance seems to have a problem with repeatability or to perform improperly, check as described below. If improper performance persists after checking, contact the local A&D dealer for repair.

Checking that the operating environment or weighing method is proper

Operating environment

- Is the weighing table solid enough?
- Is the balance level?
- Is the operating environment free from vibration and drafts?
- Is there a strong electrical or magnetic noise source such as a motor near the balance?

Weighing method

- Does the weighing pan rim touch anything? Is the weighing pan assembly installed correctly?
- Is the RE-ZERO key pressed before placing a sample on the weighing pan?
- Is the sample placed in the center of the weighing pan?
- Is the fine range breeze break ring installed for weighing with a minimum display of 0.01 mg for the GH-252 and GH-202?
- Has the balance been calibrated using the internal mass (one-touch calibration)?
- Has the balance been warmed up for one hour before weighing?

Sample and container

- Has the sample absorbed or lost moisture due to the ambient conditions such as temperature and humidity?
- Has the temperature of the container been allowed to equalize to the ambient temperature?
- Is the sample charged with static electricity?
- Is the sample of magnetic material such as iron? There are cautions about weighing magnetic materials.

Checking that the balance performs properly

- Check the balance performance using an external weight. Be sure to place the weight in the center of the weighing pan.
- Check the balance repeatability, linearity and calibrated value using external weights with a known value.

16.5. Asking for Repair

If the balance needs service or repair, contact your local A&D dealer.

The balance is a precision instrument. Use much care when handling the balance and observe the following when transporting the balance.


- Use the original packing material for transportation.
- Remove the weighing pan, pan support, breeze break ring and dust plate from the main unit.

17. Specifications

		GH-120	GH-200	GH-300	GH-202	GH-252	
Weighing capacity		120 g	220 g	320 g	220 g	250 g	
					51 g	101 g	
Maxim	num display	120.0084 g	220.0084 g	320.0084 g	220.0084 g	250.0084 g	
WIGXIII		120.000+ g	220.000+ g	020.000+ g	51.00009 g	101.00009 g	
Minimi	um display		0.1 mg		mg		
	· ·	0.1 mg			0.01 mg		
	itability	0.1	ma	0.2mg		mg	
(Stand	lard deviation)	o.r mg		0.2.119	0.02 mg	0.03 mg	
Linear	itv	±0.2	2 mg	±0.3 mg		2 mg	
					±0.03 mg	±0.10 mg	
Stabiliz	zation time	Apr	orox. 3.5 seco	onds		.5 seconds	
	(Typical at FAST)	- 4-1			Approx. 8	8 seconds	
	ivity drift,			±2 ppm/°C			
	to 30°C / 50°F to 86°F						
Operating environment		5°C to 40°C, 85%RH or less (No condensation)					
· ·	y refresh rate	5 times/second or 10 times/second					
Counting mode	Minimum unit mass	0.1 mg					
ountin	Number of						
3 -	samples	10, 25, 50 or 100 pieces					
	Minimum 100%	10.0 mg					
Percent mode	reference mass	10.0 mg					
Pe	Minimum 100% display	0.01 %, 0.1 %, 1 % (Depends on the reference mass stored.)					
Interfa	ce	RS-232C wi	th Windows C	Communicatio	n Tools Softw	are WinCT	
				300g			
Evtern	al calibration		200g	200g	20)0g	
External calibration weight		100g	100g	100g	10)0g	
		50g			5	0g	
					20g		
Weighing pan		φ90 mm					
External dimensions		217(W) x 442(D) x 316(H) mm					
Power supply &		Power consumption: Approx. 11VA (supplied to the AC adapter)					
AC adapter type		Confirm that the adapter type is correct for the local voltage and					
	1 - 91 -	power receptacle type.					
Mass		Approx. 8.2 kg					

- □ The operating environment does not include excessive change of ambient temperature, humidity, vibration, drafts, magnetic fields and static electricity.
- The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.


17.1.External Dimensions

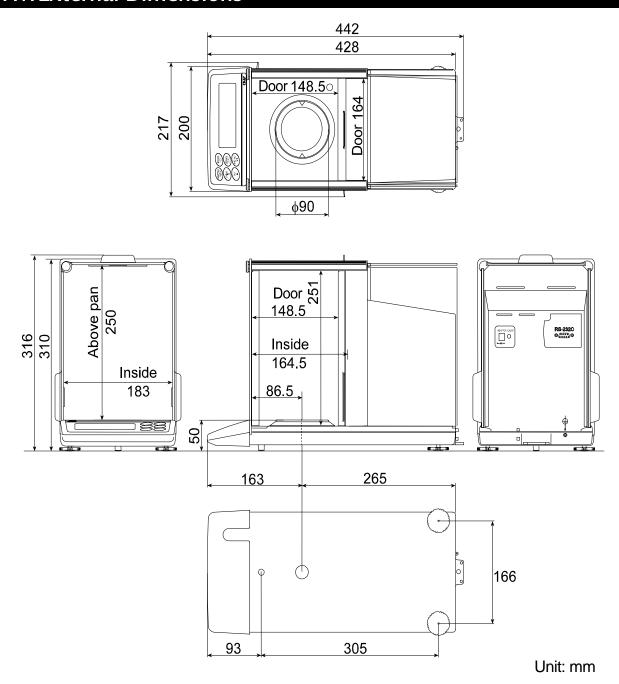

17.2. Options and Peripheral Equipment


GH-02 USB interface

- The interface to connect a balance to a personal computer, and is used to transmit the balance weight data to the personal computer via USB. Applicable OS is Windows 98 or later.
- As the Windows standard driver is used to transmit the weighing data, complicated installation of a dedicated driver is not necessary.
- The balance weighing data can be transmitted to applications such as Excel, Word and memo pad for Windows automatically.
- To perform bi-directional communications using WinCT or to output GLP data to a personal computer by using the personal computer USB interface, use the AX-USB-9P USB converter.

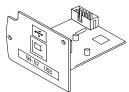

GH-08 Ethernet Interface

- The interface to connect the balance to a LAN.
- Accessory: The "WinCT-Plus" data communication software.
 - The software can acquire data from multiple balances connected to a LAN.
 - The software can control these balances with commands.
 - The software can acquire data transmitted from balances. Example: When pressing the | PRINT | key of the balance, data is output and is acquired by the computer.
 - The stored data can be used with Microsoft EXCEL (if installed).

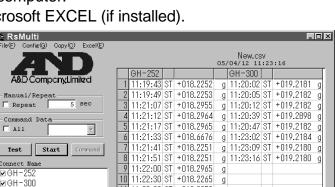


11:22:33 ST +018.2252 11:22:40 ST +018.2239


AD-8121B Printer

- Compact dot-matrix printer
- Statistical function, clock and calendar function, interval print function, graphic print function, dump print mode
- 5 x 7 dots, 16 characters per line
- Print paper (AX-PP143, 45 (W) x 50 (L) mm, ø65 mm)
- AC adapter or alkaline battery



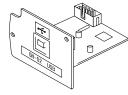

AD-1653 Density Determination Kit

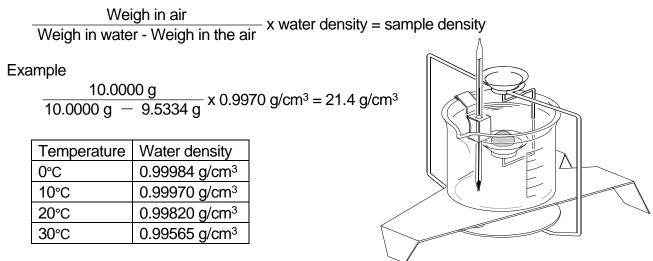


AD-1683 DC Static Eliminator

- The eliminator is able to dissipate static electricity on a charged sample.
- The eliminator does not use a ventilator and can be used in precision powder weighting.



AX-USB-9P-EX USB Converter


- Adds a COM port to a PC.
- Enables bi-directional communication between the PC and the balance when a USB driver is installed.
- Can use serial communication software such as WinCT on a PC without COM ports.
- An RS-232C cable is provided to connect the USB converter to the balance.

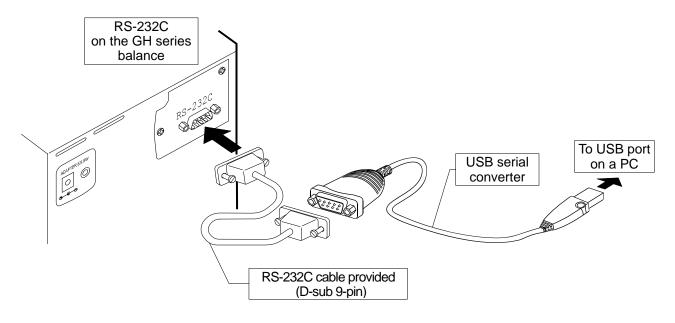

AX-T-314A-S External key Plug

This plug produces the same operations as pressing the RE-ZERO and PRINT keys. It enables remote control of the balance using an external switch.


This operation must connect the GND line to the PRINT or RE-ZERO line for at least 100 miliseconds.

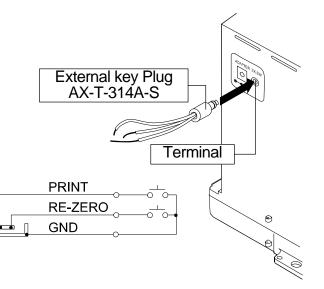

AX-SW128 Foot switch

The switch can be used to connect the external key plug and used to the RE-ZERO or PRINT keys.

Note When the foot switch is to be used, prepare the AX-SW128 and AX-T-314A-S beforehand.

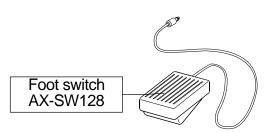

Other options

Order code	Name and description		
	Rechargeable Battery		
AD-1682	This option allows use of the balance in a place where AC power is not available.		
	Electrostatic field meter		
AD-1684	 This option measures the amount of the static charge on the sample, tare or peripheral equipment and displays the result. If those are found to be charged, discharge them using the AD-1683 DC static eliminator. 		
	Anti-vibration table		
	 This option can be used when the balance has difficulty stabilizing, due to vibration coming from the floor. 		
AD-1685	As the table has a high vibration absorbing property, key operations on the balance may make the top of the table tilt slightly, causing a weighing error. To avoid this, use the AD-8922 remote controller to control the balance.		
	Tweezers for calibration weight		
AD-1689	This option is the same as the tweezers provided as a standard accessory and is used when calibrating the balance using an external weight.		
	LAN Converter		
AD-8526	This option can be used to connect the RS-232C interface of the balance to the Ethernet (LAN) port of a computer. This allows management of the balance weighing data with a computer connected to a network.		
	Remote Display		
AD-8920	This option can be connected to the balance using the RS-232C interface.		
	Remote Controller		
AD-8922	 This option can be connected to the balance using the RS-232C interface and can control the balance remotely. 		
	 Various options such as comparator output or analog output are available. 		

18. Terms/Index

18.1.Terms

Calibration Adjustment of the balance so that it can weigh accurately.

Calibration weight A weight used for calibration

stored.

Digit The minimum weighing value available. Used for the balance, one digit

is the smallest mass that can be displayed.

Environment Ambient conditions such as vibration, drafts, temperature, static

electricity and magnetic fields which affect the weighing operation.

GLP Good Laboratory Practice.

GMP Good Manufacturing Practice

Internal mass Built-in calibration weight

ISO International Organization for Standardization

Repeatability Variation in measured values obtained when the same mass is placed

and removed repetitively. Usually expressed as a standard deviation. Example: Standard deviation = 1 digit: This means that measured values, obtained when the same sample is placed and removed

repetitively, fall within ±1 digit in the frequency of about 68%.

Re-zero To set the display to zero.

Sensitivity drift An affect that a change in temperature causes to the weighing data.

Expressed as a temperature coefficient.

Example: Temperature coefficient = 2 ppm/°C : If a load is 10 g and the temperature changes by 10°C, the value displayed changes by the

following value.

0.0002%/°C x 10°C x 10g = 0.0002 g

Stabilization time Time required after a sample being placed, until the stabilization

indicator illuminates and the weighing data is displayed.

Tare To cancel the mass value of a container which is not to be included in

the weighing data. Normally, refers to an operation of placing a

container and setting the display to zero.

Target mass An external weight used for calibration test

Zero point A weighing reference point. Usually refers to the value displayed when

nothing is on the weighing pan.

18.2. Index

%	Unit of percent mode	20	LL out	Calibration test		
\overline{lack}	0.41 1	04	CH O	Response error		71
CAL	CAL key 10,	31	CL A97	Clock		34
$\overline{\mathbb{G}}$			[Lr	Initializing the balance		30
MODE	MODE key	10	[and	Condition21, 3		
1/40			[rLF	Terminator 3		
(I/O ON:OFF	ON OFF key	10	[5] in	Internal mass correction	,	36
	·					11
(PRINT)	PRINT key 10,	31	[ES	CTS, RTS control	26	T I
	-,					
(1/10d RANGE)	RANGE key10,	31	CŢS		• • • • • •	00
771102	10 11 10 L 10 J 10,	0.	- d -	Weighing data with calender		52
→0/T← RE-ZERO	RE-ZERO key10,	21	dRER -	Data memory		
	•		Data number	er		
) \$(\$(\$)(Data number 10, 41,	53	Digit			80
▼	Interval memory mode 10,		d-no	Data number output		35
; ≮			dout	Data output34, 3		
,	Prior notice indicator10,		dS Fnc	Density function		
◀	Process indicator	10	d-E	Weighing data		52
RESPONSE FAST	MID. SLOW Response indicator	10				
	Space mark 42, 47, 48, 49, 50,		Dump print	Majahina non Error	•••••	33 70
_			- E	Weighing pan Error	•••••	70
0	Stabilization indicator 10, 16,		t	Overload error		
◀	Standby indicator	10	EC,E00	Communications error		
			EC,E01	Undefined command error		
			EC,E02	Not ready		72
A&D format		39	EC,E03	Timeout error		72
	6,		EC,E04	Excess characters error		
	on		EC,E06	Format error		
	Density determination kit		EC,E07	Parameter setting error		
			EC,E11	Stability error		
AD-1682	Rechargeable battery	79	,			
AD-1683	DC static eliminator		EC,E16	Internal mass error		
AD-1684	Electrostatic field meter		EC,E17	Internal mass error		
	Anti-vibration table	79	EC,E20	Calibration weight error		
AD-1689	Tweezers for calibration weight	79	EC,E21	Calibration weight error		
AD-8121B	Printer	76	Er[d	AK, Error code	;	36
AD-8526	LAN converter	79	Error I	Stability error		70
AD-8920	Remote display	79	Error 6	Internal mass error		70
AD-8922	Remote controller	79	Error 7	Internal mass error		70
AK code			Frror code			
	Auto print difference 34, 38,		External we	ight2	23	20 20
			FAST	2		
	Auto print polarity		FCC			
Ar-d	Zero after output		FUL			
RE-F	Auto feed			Memory full53, 5		
AX-SW128	Foot switch		GH-02	USB interface		
	S External key plug		GH-08	Ethernet Interface		
	-EX USB converter	77	GLP			
6RSFnc	Environment display31,	34	GMP			80
6PS	Baud rate		H 15	Calibration report		52
	Data bit, parity bit		ıd	ID number setting		36
[5 in	Internal mass correction		ID Number	2	41.	46
-ERL E	Calibration weight error		inFa	GLP output		
-LIIL L [AL E			int	Interval time	•••••	びと
	Calibration weight error		KF format			
	One-touch calibration			Liquid donoity input	'	4 U
[RL out	Calibration using an external weight	26	Ld in	Liquid density input	•••••	30
Calibration	23, 27,		Lo	Sample mass error		11
	est		MID.	·		21
[[in	Calibration test	25	MT format		٠ '	40

Numerical for	ormat	40	SLOW			21
Options			5Pd	Display refresh rate	34,	37
out	Report output55,	56	Stabilization	time	74,	80
P5	Internal mass correction	30	SERrE	Title block		51
PnE	Decimal point34,	37	5t-b	Stability band width	34,	37
P-on	Auto display-ON34,	37	5-Ed	Time/Date output		35
PrŁ	Data output mode 34, 38,					
PUSE	Data output pause35,	62	Target mass			80
rEcEnd	End block	51	Erc	Zero tracking	34,	37
Repeatabilit	y74,	80	Ł-UP	Timeout		36
Re-zero	•	80	£ 4PE	Data format	35,	62
RsCom		63	PCS	Unit of counting mode		18
RsKey		63	Unit	1	0, 11,	43
rtc	Clock battery error	71	טה ול	Unit		36
RTS		65	WinCT	Communication Tools		63
5- ıd	ID number output35,		Zero point			80
5 iF	Serial interface	35	-			

MEMO

	_
	_
	_
	_

MEMO

-	